

Hoechst 33342活细胞染色液(100X)

产品编号	产品名称	包装
C1028	Hoechst 33342活细胞染色液(100X)	0.5ml

产品简介:

- Hoechst 33342活细胞染色液(100X) (Hoechst 33342 Staining Solution for Live Cells, 100X), 是一种适用于活细胞细胞核染色的溶液。仅需染色10分钟就能用荧光显微镜观察到非常明亮的细胞核蓝色荧光染色。
- 本产品也适用于固定后的细胞或组织的细胞核染色。
- Hoechst 33342, 也称bisBenzimide H 33342或HOE 33342。分子式为 $C_{27}H_{28}N_6O \cdot 3HCl \cdot 3H_2O$, 分子量为615.99, CAS Number 23491-52-3。Hoechst 33342的最大激发波长为346nm, 最大发射波长为460nm; Hoechst 33342和双链DNA结合后, 最大激发波长为350nm, 最大发射波长为461nm。
- Hoechst 33342是一种可以穿透细胞膜的蓝色荧光染料, 对细胞的毒性较低。
- Hoechst 33342染色常用于细胞凋亡检测, 染色后用荧光显微镜观察或流式细胞仪检测。Hoechst 33342也常用于普通的细胞核染色, 或常规的DNA染色。
- 本产品用于活细胞染色的效果图参考图1。

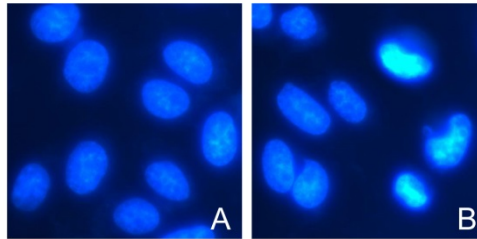


图1. 正常培养的HeLa细胞用本产品染色后的实拍效果图。图A为正常细胞Hoechst 33342的染色效果图。图B中除了正常细胞外, 还清晰可见呈现致密浓染的凋亡细胞。

包装清单:

产品编号	产品名称	包装
C1028	Hoechst 33342活细胞染色液(100X)	0.5ml
—	说明书	1份

保存条件:

-20°C避光保存, 一年有效。

注意事项:

- 本产品浓度经过碧云天的优化, 确保可以满足活细胞染色及各种常规染色的需要。如需使用特定浓度的Hoechst 33342, 请选购碧云天的粉末装Hoechst 33342 (C1022)。
- 荧光染料都存在淬灭的问题, 建议染色后尽量当天完成检测。
- 为减缓荧光淬灭可以使用抗荧光淬灭封片液。抗荧光淬灭封片液(P0126)可以向碧云天订购。
- Hoechst 33342对人体有害, 操作时请小心, 并注意有效防护以避免直接接触人体或吸入体内。
- 本产品仅限于专业人员的科学研究用, 不得用于临床诊断或治疗, 不得用于食品或药品, 不得存放于普通住宅内。
- 为了您的安全和健康, 请穿实验服并戴一次性手套操作。

使用说明:

1. 对于活细胞或培养的组织:

- 对于培养器皿内的活细胞或组织, 在培养液中均匀滴加适量的Hoechst 33342活细胞染色液(100X)至终浓度为1X, 轻柔混匀。例如对于十二孔板中培养的细胞, 每孔有1ml的培养液, 每孔需要加入10 μ l Hoechst 33342活细胞染色液(100X)。培养液中的血清、酚红等都不会对染色产生干扰。
- 在适宜于细胞培养的温度孵育10分钟。吸除含染料的培养液, 用培养液或PBS洗涤2-3次即可在荧光显微镜下观察。细胞发生凋亡时, 会看到凋亡细胞的细胞核呈致密浓染, 或呈碎块状致密浓染。具体示例见图1。如果对于染色效果的要求不是特别高, 染色5分钟, 用培养液或PBS洗涤1次即可在荧光显微镜下观察。为避免凋亡细胞随培养液或PBS被吸除, 在吸除液体前, 对于用多孔板中培养的细胞, 最好用多孔板离心机离心一下以充分沉淀那些已经漂浮起来的凋亡细胞。

2. 对于固定的细胞或组织:

- a. 对于细胞或组织样品, 固定后, 适当洗涤去除固定剂。随后如果需要免疫荧光染色, 则先进行免疫荧光染色, 染色完后再按后续步骤进行Hoechst 33342染色。如果不需要进行其它染色, 则直接进行后续的Hoechst 33342染色。
- b. 对于贴壁细胞或组织切片, 加入少量用免疫染色洗涤液、PBS或生理盐水稀释至1X的本产品, 覆盖住样品即可。对于悬浮细胞, 至少加入待染色样品体积3倍的染色液, 混匀。室温放置3-5分钟。
- c. 吸除Hoechst 33342染色液, 用免疫染色洗涤液、PBS或生理盐水洗涤2-3次, 每次3-5分钟。
- d. 直接在荧光显微镜下观察或封片后荧光显微镜下观察。细胞发生凋亡时, 会看到凋亡细胞的细胞核呈致密浓染, 或呈碎块状致密浓染。

相关产品:

产品编号	产品名称	包装
C1002	DAPI	5mg/ml×0.2ml
C1005	DAPI 染色液	10ml
C1006	DAPI 染色液	50ml
C1011	Hoechst 33258	10mg
C1017	Hoechst 33258 染色液	10ml
C1018	Hoechst 33258 染色液	50ml
C1022	Hoechst 33342	10mg
C1025	Hoechst 33342 染色液	10ml
C1026	Hoechst 33342 染色液	50ml
C1027	Hoechst 33342 活细胞染色液(100X)	0.1ml
C1028	Hoechst 33342 活细胞染色液(100X)	0.5ml
C1029	Hoechst 33342 活细胞染色液(100X)	3ml

使用本产品的文献:

1. You F, Li Q, Jin G, Zheng Y, Chen J, Yang H.Genistein protects against A β 25-35 induced apoptosis of PC12cells through JNK signaling and modulation of Bcl-2 family messengers.BMC Neurosci . 2017 Jan 12;18(1):12
2. Wang Z, Liao K, Zuo W, Liu X, Qiu Z, Gong Z, Liu C, Zeng Q, Qian Y, Jiang L, Bu Y, Hong S, Hu G.Depletion of NFB1/MDC1 Induces Apoptosis in Nasopharyngeal Carcinoma Cells Through the p53-ROS-Mitochondrial Pathway.Oncol Res . 2017 Jan 2;25(1):123-136
3. Hu Y, Xiong LL, Zhang P, Wang TH.Microarray expression profiles of genes in lung tissues of rats subjected to focal cerebral ischemia-induced lung injury following bone marrow-derived mesenchymal stem cell transplantation.Int J Mol Med . 2017 Jan;39(1):57-70
4. Zhang Y, Wu JY, Weng LH, Li XX, Yu LJ, Xu Y.Valproic acid protects against MPP+-mediated neurotoxicity in SH-SY5Y Cells through autophagy.Neurosci Lett . 2017 Jan 18;638:60-68
5. Fan Y, Liu J, Liu D, Zhou Z, Bao Y, Wang J, Zhao Q, Xu Y.NSCA-1-a novel N-substituted coumalamide derivative-increases Adriamycin sensitivity in HepG2/adriamycin cells through modulating Akt/GSK-3 β signaling and p53-dependant apoptotic pathway.ENVIRON TOXICOL CHEM . 2017 Jan;49:1-7
6. Chang S, Ruan WC, Xu YZ, Wang YJ, Pang J, Zhang LY, Liao H, Pang T.The natural product 4,10-aromadendranediol induces neuritogenesis in neuronal cells in vitro through activation of the ERK pathway.Acta Pharmacol Sin . 2017 Jan;38(1):29-40
7. Wang Y1, Qiu Y1, Yin S1, Zhang L1, Shi K1, Gao H1, Zhang Z1, He Q1.A functional nanocarrier that copenetrates extracellular matrix and multiple layers of tumor cells for sequential and deep tumor autophagy inhibitor and chemotherapeutic delivery.Autophagy . 2017 Feb;13(2):359-370
8. Wen Y, He W, Jiang M, Zeng M, Cai L.Deriving cells expressing markers of female germ cells from premature ovarian failure patient-specific induced pluripotent stem cells.Regen Med . 2017 Mar;12(2):143-152
9. Zhai Y1, Li Y2, Wang Y3, Cui J2, Feng K4, Kong X2, Chen L5.Psoralidin, a prenylated coumestan, as a novel anti-osteoporosis candidate to enhance boneformation of osteoblasts and decrease bone resorption of osteoclasts.Eur J Pharmacol . 2017 Apr 15;801:62-71
10. Wu J, Yang X, Lu H, Liu L, Xu B, Zheng S, Yu B, Jie K, Wan F.Inhibitory effect and mechanism of exogenous mammalian sterile 20-like kinase 1 on the growth of human colorectal cancer.Oncol Lett . 2017 Apr;13(4):2656-2664
11. Guo Y, Song Z, Zhou M, Yang Y, Zhao Y, Liu B, Zhang X.Infiltrating macrophages in diabetic nephropathy promote podocytes apoptosis via TNF- α -ROS-p38MAPK pathway.ONCOTARGET . 2017 Jun 7;8(32):53276-53287
12. Sun L, Li Y, Jia X, Wang Q, Li Y, Hu M, Tian L, Yang J, Xing W, Zhang W, Wang J, Xu H, Wang L, Zhang D, Ren H.Neuroprotection byIFN- γ via astrocyte-secreted IL-6 in acute neuroinflammation.ONCOTARGET . 2017 Jun 20;8(25):40065-40078
13. Xiong LL, Liu F, Lu BT, Zhao WL, Dong XJ, Liu J, Zhang RP, Zhang P, Wang TH.Bone Marrow Mesenchymal Stem-Cell Transplantation Promotes Functional ImprovementAssociated with CNTF-STAT3 Activation after Hemi-Sectioned Spinal Cord Injury in Tree Shrews.Front Cell Neurosci . 2017 Jun 28;11:172
14. Wu X, Gao H, Sun W, Yu J, Hu H, Xu Q, Chen X.Nepetoidin B,a Natural Product, Inhibits LPS-stimulated Nitric Oxide Production via Modulation of iNOS Mediated by NF- κ B/MKP-5 Pathways.Phytother Res . 2017 Jul;31(7):1072-1077
15. Yan X, Xun M, Dou X, Wu L, Han Y, Zheng J.Regulation of Na $^{+}$ -K $^{+}$ -ATPase effected high glucose-induced myocardial cell injury through c-Srcdependent NADPH oxidase/ROS pathway.Exp Cell Res . 2017 Aug 15;357(2):243-251
16. Liu JC, Lai FN, Li L, Sun XF, Cheng SF, Ge W, Wang YF, Li L, Zhang XF, De Felici M, Dyce PW, Shen W.Di (2-ethylhexyl) phthalate exposure impairs meiotic progression and DNA damage repair in fetal mouse oocytes in vitro.Cell Death Dis . 2017 Aug 3;8(8):e2966.
17. Liu XY, Zhang LJ, Chen Z, Liu LB.The PTEN inhibitor bpV(pic) promotes neuroprotection against amyloid β -peptide(25-35)-induced oxidative stress and neurotoxicity.Neurol Res . 2017 Aug;39(8):758-765
18. Shan ZL, Zhong L, Xiao CL, Gan LG, Xu T, Song H, Yang R, Li L, Liu BZ.Shikonin suppresses proliferation and induces apoptosis in human leukemia NB4

cells through modulation of MAPKs and c Myc. *Mol Med Rep* . 2017 Sep;16(3):3055-3060

19. Hao J, Yang Z, Wang L, Zhang Y, Shu Y, Jiang L, Hu Y, Lv W, Dong P, Liu Y. Downregulation of BRD4 inhibits gallbladder cancer proliferation and metastasis and induces apoptosis via PI3K/AKT pathway. *Int J Oncol* . 2017 Sep;51(3):823-831. (IF 3.899)
20. Liu X, Qiu Z, Wang Z, Zuo W, Gong Z, Liu C, Zeng Q, Qian Y, Jiang L, Li Y, Bu Y, Hu G. NFB1/MDC1 participates in the regulation of proliferation and apoptosis in human laryngeal squamous cell carcinoma. *Clin Transl Oncol* . 2017 Sep 18
21. Ying C, Xiao BD, Qin Y, Wang BR, Liu XY, Wang RW, Fang L, Yan H, Zhou XM, Wang YG. GOLPH2-regulated oncolytic adenovirus, GD55, exerts strong killing effect on human prostate cancer stem-like cells in vitro and in vivo. *Acta Pharmacol Sin* . 2017 Sep 7
22. Liang J, Wang WF, Xie S, Zhang XL, Qi WF, Zhou XP, Hu JX, Shi Q, Yu RT. β -transducin repeat-containing E3 ubiquitin protein ligase inhibits migration, invasion and proliferation of glioma cells. *Oncol Lett* . 2017 Sep;14(3):3131-3135
23. Geng YD, Zhang C, Lei JL, Yu P, Xia YZ, Zhang H, Yang L, Kong LY. Walsuronoid B induces mitochondrial and lysosomal dysfunction leading to apoptotic rather than autophagic cell death via ROS/p53 signaling pathways in liver cancer. *Biochem Pharmacol* . 2017 Oct 15;142:71-86
24. Chen GH, Xu CS, Zhang J, Li Q, Cui HH, Li XD, Chang LP, Tang RJ, Xu JY, Tian XQ, Huang PS, Xu J, Jin C, Yang YJ. Inhibition of miR-128-3p by Tongxinluo Protects Human Cardiomyocytes from Ischemia/reperfusion Injury via Upregulation of p70s6k1/p-p70s6k1. *Front Pharmacol* . 2017 Oct 30;8:775
25. Liu W, Wang X, Liu Z, Wang Y, Yin B, Yu P, Duan X, Liao Z, Chen Y, Liu C, Li X, Dai Y, Tao Z. SGK1 inhibition induces autophagy-dependent apoptosis via the mTOR-Foxo3a pathway. *BRIT J CANCER* . 2017 Oct 10;117(8):1139-1153
26. Li Y, Sun Y, Cao T, Su Q, Li Z, Huang M, Ouyang R, Chang H, Zhang S, Miao YA. A cation-exchange controlled core-shell MnS@Bi2S3 theranostic platform for multimodal imaging guided radiation therapy with hyperthermia boost. *Nanoscale* . 2017 Oct 5;9(38):14364-14375
27. Qian L, Li X, Ye P, Wang G, Dai W, Liu Y, Gao Q, Shen G. Oxymatrine induces apoptosis and inhibits invasion in Gallbladder carcinoma via PTEN/PI3K/AKT pathway. *Cytotechnology* . 2017 Nov 23
28. Zhao C, Yao X, Chen X, Wu W, Xi F, Yang G, Yu T. Knockdown of ubiquitin D inhibits adipogenesis during the differentiation of porcine intramuscular and subcutaneous preadipocytes. *CELL PROLIFERAT* . 2017 Nov 23
29. Yang LL, Zhao Y1, Luo SM1, Ma JY1, Ge ZJ1, Shen W1, Yin S2. Toxic Effects and Possible Mechanisms of Hydrogen Sulfide and/or Ammonia on Porcine Oocyte Maturation in vitro. *Toxicol Lett* . 2017 Dec 29. pii: S0378-4274(17)31529-1
30. Xiao B, Qin Y, Ying C, Ma B, Wang B, Long F, Wang R, Fang L, Wang Y. Combination of oncolytic adenovirus and luteolin exerts synergistic antitumor effects in colorectal cancer cells and a mouse model. *Mol Med Rep* . 2017 Dec;16(6):9375-9382
31. Li X, Yan X, Wang Y, Wang J, Zhou F, Wang H, Xie W, Kong H. NLRP3 inflammasome inhibition attenuates silica-induced epithelial to mesenchymal transition (EMT) in human bronchial epithelial cells. *Exp Cell Res* . 2017 Dec 16. pii: S0014-4827(17)30668-7
32. Xu Z, Zheng M, Zhang L, Zhang X, Zhang Y, Cui X, Gong X, Xi R, Bai R. Effect of ATP and Bax on the apoptosis of Eimeria tenella host cells. *BMC Vet Res* . 2017 Dec 28;13(1):399
33. Sun Y, Wang C, Meng Q, Liu Z, Huo X, Sun P, Sun H, Ma X, Peng J, Liu K. Targeting P-glycoprotein and SORCIN: Dihydropyridinyl strengthens anti-proliferative efficiency of adriamycin via MAPK/ERK and Ca²⁺-mediated apoptosis pathways in MCF-7/ADR and K562/ADR. *J Cell Physiol* . 2018 Apr;233(4):3066-3079
34. Guo Q, Zheng M, Xu Y, Wang N, Zhao W. miR-384 induces apoptosis and autophagy of non-small cell lung cancer cells through the negative regulation of Collagen α -1(X) chain gene. *BIOSCIENCE REP* . 2019 Feb 1;39(2). pii: BSR20181523
35. Ayman M Mahmoud, Fiona L Wilkinson, Mansur A Sandhu, Julia M Dos Santos, M Yvonne Alexander. Modulating Oxidative Stress in Drug-Induced Injury and Metabolic Disorders: The Role of Natural and Synthetic Antioxidants. *Oxid Med Cell Longev* . 2019 Jun 27;2019:3206401.;doi: 10.1155/2019/3206401
36. Zhang J, Si H, Li B, Zhou X, Zhang J. Myrislignan Exhibits Activities Against Toxoplasma gondii RH Strain by Triggering Mitochondrial Dysfunction. *Front Microbiol* . 2019 Sep 18;10:2152
37. Tang LX, Su SF, Wan Q, He P, Xiang Y, Cheng XM. Novel long non-coding RNA LBX2-AS1 indicates poor prognosis and promotes cell proliferation and metastasis through Notch signaling in non-small cell lung cancer. *EUR REV MED PHARMACO* . 2019 Sep;23(17):7419-7429
38. Jiyue Ge, Likai Mao, Weiqiang Xu, Wenge Fang, Ningning Wang, Dawen Ye, Zhuang Dong, Han Guan, Chao Guan. miR-103a-3p Suppresses Cell Proliferation and Invasion by Targeting Tumor Protein D52 in Prostate Cancer. *J Invest Surg* . 2020 Mar 13;1-9.;doi: 10.1080/08941939.2020.1738602
39. Lingqi Zhou, Hai Tang, Fang Wang, Shanshan Ou, Tong Wu, Yinchao Fang, Jie Xu, Kaihua Guo. Cyclovirobuxine D inhibits cell proliferation and migration and induces apoptosis in human glioblastoma multiforme and low-grade glioma. *Oncol Rep* . 2020 Mar;43(3):807-816.;doi: 10.3892/or.2020.7459
40. Peng Chen, Xuejie Li, Ruonan Zhang, Shuiping Liu, Yu Xiang, Mingming Zhang, Xiaying Chen, Ting Pan, Lili Yan, Jiao Feng, Ting Duan, Da Wang, Bi Chen, Ting Jin, Wengang Wang, Liuxi Chen, Xingxing Huang, Wenzheng Zhang, Yitian Sun, Guohua Li, Lingpan Kong, Xiaohui Chen, Yongqiang Li, Zuyi Yang, Qin Zhang, Lvjia Zhuo, Xinbing Sui, Tian Xie. Combinative treatment of β -elemene and cetuximab is sensitive to KRAS mutant colorectal cancer cells by inducing ferroptosis and inhibiting epithelial-mesenchymal transformation. *Theranostics* . 2020 Apr 6;10(11):5107-5119.;doi: 10.7150/thno.44705
41. Lei Lai, Yizhou Xu, Lan Kang, Jianmin Yang, Gangjie Zhu. LncRNA KCNQ1OT1 contributes to cardiomyocyte apoptosis by targeting FUS in heart failure. *Exp Mol Pathol* . 2020 Aug;115:104480.;doi: 10.1016/j.yexmp.2020.104480
42. Li Mao, Mei-Ling Zuo, Ai-Ping Wang, Ying Tian, Li-Chen Dong, Tao-Ming Li, Da-Bin Kuang, Gui-Lin Song, Zhong-Bao Yang. Low expression of miR-532-3p contributes to cerebral ischemia/reperfusion oxidative stress injury by directly targeting NOX2. *Mol Med Rep* . 2020 Sep;22(3):2415-2423.;doi: 10.3892/mmr.2020.11325
43. Runqing Geng, Yuanyuan Ren, Rong Rao, Xi Tan, Hong Zhou, Xiangliang Yang, Wei Liu, Qunwei Lu. Titanium Dioxide Nanoparticles Induced HeLa Cell Necrosis under UVA Radiation through the ROS-mPTP Pathway. *Nanomaterials (Basel)* . 2020 Oct 15;10(10):2029.;doi: 10.3390/nano10102029
44. Ying Hua, Jiali Wu, Muqing Fu, Jinyue Liu, Xiaoxia Li, Bao Zhang, Wei Zhao, Chengsong Wan. Enterohemorrhagic Escherichia coli Effector Protein EspF Interacts With Host Protein ANXA6 and Triggers Myosin Light Chain Kinase (MLCK)-Dependent Tight Junction Dysregulation. *Front Cell Dev Biol* . 2020 Dec 23;8:613061.;doi: 10.3389/fcell.2020.613061
45. Gu Y, Cai R, Zhang C, Xue Y, Pan Y, Wang J, Zhang Z. miR-132-3p boosts caveolae-mediated transcellular transport in glioma endothelial cells by targeting PTEN/PI3K/PKB/Src/Cav-1 signaling pathway. *ANN ONCOL*

Version 2021.09.01