



碧云天网站



微信公众号

碧云天生物技术/Beyotime Biotechnology

订货热线: 400-168-3301或800-8283301

订货e-mail: order@beyotime.com

技术咨询: info@beyotime.com

网址: http://www.beyotime.com

化学发光法 EMSA 试剂盒

产品编号	产品名称	包装
GS009	化学发光法EMSA试剂盒	100次

产品简介:

- 化学发光法EMSA试剂盒(Chemiluminescent EMSA Kit)是一种通过Streptavidin-HRP及后续的BeyoECL Moon试剂来实现化学发光检测Biotin标记的EMSA探针的检测试剂盒。同时本试剂盒也提供了EMSA检测所需的结合缓冲液和上样缓冲液, 及一些关键的相关试剂, 可以实现非同位素的EMSA检测。
- 本试剂盒采用了高质量的Streptavidin-HRP Conjugate, HRP和Streptavidin共价交联的比例大于3, 这样比采用Streptavidin和Biotin-HRP conjugate两种试剂进行检测更方便, 并且灵敏度更高。
- 本试剂盒采用了非特异性结合比avidin更低的streptavidin, 使检测结果背景更低灵敏度更高。
- 本试剂盒和碧云天的各种生物素标记EMSA探针可以配套使用。
- 本试剂盒没有提供生物素探针标记相关的试剂, 其它特定的生物素标记EMSA探针的制备可以使用碧云天的EMSA探针生物素标记试剂盒(GS008)。
- EMSA/Gel-Shift 结合缓冲液(5X)中含有poly(dI-dC)等有效成分。其中poly(dI-dC)的浓度经过优化, 可以很好的消除蛋白和标记探针间的非特异性结合, 同时又不会减弱目的转录因子和标记探针间的结合。
- 本试剂盒可以用于100个蛋白和探针的结合反应, 并足够检测至少10块有生物素标记EMSA探针的膜。

包装清单:

产品编号	产品名称	包装
GS009-1	EMSA/Gel-Shift结合缓冲液(5X)	200μl
GS009-2	EMSA/Gel-Shift上样缓冲液(无色, 10X)	200μl
GS009-3	EMSA/Gel-Shift上样缓冲液(蓝色, 10X)	200μl
GS009-4	BeyoECL Moon A液	55ml
GS009-5	BeyoECL Moon B液	55ml
GS009-6	Streptavidin-HRP Conjugate	100μl
GS009-7	封闭液	380ml
GS009-8	洗涤液(5X)	250ml
GS009-9	检测平衡液	250ml
—	说明书	1份

保存条件:

GS009-1至GS009-3和GS009-6在-20℃保存, 其余可4℃保存。如果长期不用, 整个试剂盒可-20℃保存, -20℃可以保存更长时间。

注意事项:

- 需自备带正电荷尼龙膜, 以及凝胶电泳时所需的相关试剂。带正电荷尼龙膜(FFN10/FFN11/FFN13/FFN15)可以向碧云天订购。
- 如果需要使用更多的封闭液或洗涤液, 可另外单独订购GS009B封闭液或GS009W洗涤液(5X)。
- BeyoECL Moon A液和B液均对人体有害, 操作时请小心, 并注意有效防护以避免直接接触人体或吸入体内。
- 本产品仅限于专业人员的科学研究用, 不得用于临床诊断或治疗, 不得用于食品或药品, 不得存放于普通住宅内。
- 为了您的安全和健康, 请穿实验服并戴一次性手套操作。

使用说明:

1. 探针的标记:

可以直接选购碧云天的生物素标记EMSA探针, 或使用碧云天的EMSA探针生物素标记试剂盒(GS008)或其它合适的试剂盒进行EMSA探针的生物素标记。

2. 探针的纯化和检测:

碧云天的各种生物素标记EMSA探针都已经过纯化, 可以直接使用; 对于使用碧云天的EMSA探针生物素标记试剂盒(GS008)等试剂盒标记的EMSA探针, 通常为实验简便起见, 可以不必纯化标记好的探针。有些时候, 纯化后的探针会改善

EMSA的电泳结果。详细的探针纯化和探针的标记效率的检测请参考EMSA探针生物素标记试剂盒的相关说明。

3. EMSA胶的配制:

- 准备好倒胶的模具。可以使用常规的制备蛋白电泳胶的模具(例如BioRad的常规用于蛋白电泳的制胶装置), 或其它适当的模具。最好选择可以灌制较薄胶的模具, 以便于干胶等后续操作。为得到更好的结果, 可以选择可灌制较大EMSA胶的模具。制胶前必须把制胶模具冲洗干净, 需特别注意不能有SDS残留。
- 按照如下配方配制20ml 4%的聚丙烯酰胺凝胶(注意: 使用29:1等不同比例的Acr/Bis对结果影响不大)。

TBE buffer (10X)	1.0ml
重蒸水	16.2ml
39:1 acrylamide/bisacrylamide (40%,w/v)	2ml
80% 甘油	625 μ l
10% 过硫酸铵(ammonium persulfate)	150 μ l
TEMED	10 μ l

- 按照上述顺序依次加入各种试剂, 加入TEMED前先混匀, 加入TEMED后立即混匀, 并马上加入到制胶的模具中。避免产生气泡, 并加上梳齿。如果发现非常容易形成气泡, 可以把一块制胶的玻璃板进行硅烷化处理。

4. EMSA结合反应:

- 如下设置EMSA结合反应:

阴性对照反应:

Nuclease-Free Water	7 μ l
EMSA/Gel-Shift 结合缓冲液(5X)	2 μ l
细胞核蛋白或纯化的转录因子	0 μ l
标记好的探针	1 μ l
总体积	10 μ l

探针冷竞争反应:

Nuclease-Free Water	4 μ l
EMSA/Gel-Shift 结合缓冲液(5X)	2 μ l
细胞核蛋白或纯化的转录因子	2 μ l
未标记的探针	1 μ l
标记好的探针	1 μ l
总体积	10 μ l

Super-shift反应:

Nuclease-Free Water	4 μ l
EMSA/Gel-Shift 结合缓冲液(5X)	2 μ l
细胞核蛋白或纯化的转录因子	2 μ l
目的蛋白特异抗体	1 μ l
标记好的探针	1 μ l
总体积	10 μ l

样品反应:

Nuclease-Free Water	5 μ l
EMSA/Gel-Shift 结合缓冲液(5X)	2 μ l
细胞核蛋白或纯化的转录因子	2 μ l
标记好的探针	1 μ l
总体积	10 μ l

突变探针的冷竞争反应:

Nuclease-Free Water	4 μ l
EMSA/Gel-Shift 结合缓冲液(5X)	2 μ l
细胞核蛋白或纯化的转录因子	2 μ l
未标记的突变探针	1 μ l
标记好的探针	1 μ l
总体积	10 μ l

- 按照上述顺序依次加入各种试剂, 在加入标记好的探针前先混匀, 并且室温(20-25 $^{\circ}$ C)放置10分钟, 从而消除可能发生的探针和蛋白的非特异性结合, 或者让冷探针优先反应。然后加入标记好的探针, 混匀, 室温(20-25 $^{\circ}$ C)放置20分钟。
- 加入1 μ l EMSA/Gel-Shift上样缓冲液(无色, 10X), 混匀后立即上样。注意: 有些时候溴酚蓝会影响蛋白和DNA的结合, 建议尽量使用无色的EMSA/Gel-Shift上样缓冲液。如果对于使用无色上样缓冲液在上样时感觉到无法上样, 可以在无色上样缓冲液里面添加极少量的蓝色的上样缓冲液, 至可以观察到蓝颜色即可。

5. 电泳:

- 用0.5XTBE作为电泳液。按照10V/厘米的电压预电泳10分钟。预电泳的时候如果有空余的上样孔, 可以加入少量稀释好的1X的EMSA上样缓冲液(蓝色), 以观察电压是否正常进行。
- 把混合了上样缓冲液的样品加入到上样孔内。在多余的某个上样孔内加入10 μ l稀释好的1X的EMSA/Gel-Shift上样缓冲液(蓝色), 用于观察电泳进行的情况。
- 按照10V/厘米的电压电泳。确保胶的温度不超过30 $^{\circ}$ C, 如果温度升高, 需要适当降低电压。电泳至EMSA/Gel-Shift上样缓冲液中的蓝色染料溴酚蓝至胶的下缘1/4处, 停止电泳。

6. 转膜:

- 取一和EMSA胶大小相近或略大的尼龙膜, 剪角做好标记, 用0.5XTBE浸泡至少10分钟。尼龙膜自始至终仅能使用镊子夹取, 并且仅可夹取不可能接触样品的边角处。
- 取两片和尼龙膜大小相近或略大的滤纸, 用0.5XTBE浸湿。
- 把浸泡过的尼龙膜放置在一片浸湿的滤纸上, 注意避免尼龙膜和滤纸间产生气泡。
- 非常小心地取出EMSA胶放置到尼龙膜上, 注意确保胶和膜之间没有气泡。
- 再把另外一片浸湿的滤纸放置到EMSA胶上, 注意确保滤纸和胶之间没有气泡。

- f. 采用Western时所使用的湿法电转膜装置或其它类似的电转膜装置,以0.5XTBE为转膜液,把EMSA胶上的探针、蛋白以及探针和蛋白的复合物等转移到尼龙膜上。对于大小约为10x8x0.1cm的EMSA胶,用BioRad的常用的Western转膜装置,电转时可以设置为380mA(约100V)转膜30-60分钟。如果胶较厚,则需适当延长转膜时间。转膜时需保持转膜液的温度较低,通常可以把电转槽置于4°C冷库或置于冰浴或冰水浴中进行电转,这样可以确保低温。具体的电转膜方法请参考电转膜装置的使用说明。
- g. 转膜完毕后,小心取出尼龙膜,样品面向上,放置在一干燥的滤纸上,轻轻吸掉下表面明显的液体。立即进入下一步的交联步骤,不可使膜干掉。
- 7. 交联:**
- a. 用紫外交联仪(UV-light cross-linker)选择254nm紫外波长,120mJ/cm²,交联45-60秒。如果没有紫外交联仪可以使用普通的手提式紫外灯(例如碧云天的手提紫外检测仪(EUV002)),距离膜5-10厘米左右照射3-10分钟。也可以使用超净工作台内的紫外灯,距离膜5-10厘米左右照射3-15分钟。最佳的交联时间可以使用标准品自行摸索。
- b. 交联完毕后,可以直接进入下一步检测;也可以用保鲜膜包裹后在室温干燥处存放3-5天,然后再进入下一步检测。
- c. 如果检测结果发现交联效果不佳,甚至连free probe的条带都非常微弱,可以考虑在膜干燥后参考步骤A的条件再交联一次,以进一步改善交联效果。
- 8. 化学发光法检测生物素标记的探针:**
- a. 37-50°C水浴溶解封闭液和洗涤液。
注意: 封闭液和洗涤液必须完全溶解后方可使用,封闭液和洗涤液可以在室温至50°C之间使用,但必须确保这两种溶液中均无沉淀产生,在冬天需特别注意。
- b. 取一合适的容器加入15ml封闭液,再放入交联过的含有样品的尼龙膜。在侧摆摇床或水平摇床上缓慢摇动15分钟。
- c. 取7.5μl Streptavidin-HRP Conjugate加入到15ml封闭液中(1:2000稀释),混匀备用。
- d. 去除用于尼龙膜封闭的封闭液,加入上一步中配制的15ml含有Streptavidin-HRP Conjugate的封闭液。在侧摆摇床或水平摇床上缓慢摇动15分钟。
- e. 取25ml洗涤液(5X),加入100ml重蒸水或Milli-Q级纯水,混匀配制125ml洗涤液。
- f. 将尼龙膜转移至另一装有15-20ml洗涤液的容器内,漂洗1分钟。
- g. 去除洗涤液,加入15-20ml洗涤液,在侧摆摇床或水平摇床上缓慢上洗涤5分钟。
- h. 重复步骤G三次(共洗涤四次),每次洗涤时间都约为5分钟。
- i. 将尼龙膜转移至另一装有20-25ml检测平衡液的容器内,在侧摆摇床或水平摇床上缓慢摇动5分钟。
- j. 取5ml BeyoECL Moon A液和5ml BeyoECL Moon B液混匀,配制BeyoECL Moon工作液。注意: BeyoECL Moon工作液必须现配现用。说明: 从本步骤起操作方法和注意事项同Western实验的荧光检测。
- k. 取出尼龙膜,用吸水纸吸去过多液体。立即将膜的样品面向上,放置到处于水平桌面上的洁净容器内或保鲜膜上。
- l. 在尼龙膜的表面小心加上步骤I配制好的共10ml BeyoECL Moon工作液,使工作液完全覆盖尼龙膜。室温放置2-3分钟。
- m. 取出尼龙膜,用吸水纸吸去过多液体。将尼龙膜放在两片保鲜膜或其它适当的透光薄膜中间,并固定于压片暗盒(也称片夹)内。
- n. 用X光片压片1-5分钟。可以先压片1分钟,立即显影定影,然后根据结果再调整压片时间;也可以直接分别压片30秒、1、3、5分钟或更长时间,然后一起显影定影观察结果。

使用本产品的文献:

- Wang X, Li C, Chen Y, Hao Y, Zhou W, Chen C, Yu Z. Hypoxia enhances CXCR4 expression favoring microglia migration via HIF-1 α activation. *Biochem Biophys Res Commun*. 2008 Jun 27;371(2):283-8.
- Zhang L, Chen W, Li X. A novel anticancer effect of butein: inhibition of invasion through the ERK1/2 and NF-kappa B signaling pathways in bladder cancer cells. *FEBS Lett*. 2008 Jun 11;582(13):1821-8.
- Wei Y, Weng D, Li F, Zou X, Young DO, Ji J, Shen P. Involvement of JNK regulation in oxidative stress-mediated murine liver injury by microcystin-LR. *Apoptosis*. 2008 Aug;13(8):1031-42.
- Mao XM, Zhou Z, Cheng LY, Hou XP, Guan WJ, Li YQ. Involvement of SigT and RstA in the differentiation of *Streptomyces coelicolor*. *FEBS Lett*. 2009 Oct 6;583(19):3145-50.
- Lu J, Zhao J, Liu K, Zhao J, Yang H, Huang Y, Qin Z, Bai R, Li P, Ma J, Yan W, Zhao M, Dong Z. MAPK/ERK1/2 signaling mediates endothelial-like differentiation of immature DCs in the microenvironment of esophageal squamous cell carcinoma. *Cell Mol Life Sci*. 2010 Jun;67(12):2091-106.
- Ding RY, Zhao DM, Zhang ZD, Guo RX, Ma XC. Pretreatment of Rho kinase inhibitor inhibits systemic inflammation and prevents endotoxin-induced acute lung injury in mice. *J Surg Res*. 2011 Dec;171(2):e209-14.
- Xu J, Wang B, Xu Y, Sun L, Tian W, Shukla D, Barod R, Grillari J, Grillari-Voglauer R, Maxwell PH, Esteban MA. Epigenetic regulation of HIF-1 α in renal cancer cells involves HIF-1 α /2 α binding to a reverse hypoxia-response element. *Oncogene*. 2012 Feb 23;31(8):1065-72.
- Gu L, Liang X, Wang L, Yan Y, Ni Z, Dai H, Gao J, Mou S, Wang Q, Chen X, Wang L, Qian J. Functional metabotropic glutamate receptors 1 and 5 are expressed in murine podocytes. *Kidney Int*. 2012 Mar;81(5):458-68.
- Peng X, Wang Y, Kolli S, Deng J, Li L, Wang Z, Raj JU, Gou D. Physical and functional interaction between the ID1 and p65 for activation of NF κ B. *Am J Physiol Cell Physiol*. 2012 Aug 1;303(3):C267-77.
- Liu SS, Zheng HX, Jiang HD, He J, Yu Y, Qu YP, Yue L, Zhang Y, Li Y. Identification and characterization of a novel gene, c1orf109, encoding a CK2 substrate that is involved in cancer cell proliferation. *J Biomed Sci*. 2012 May 1;19(1):49.
- Li F, Li C, Zhang H, Lu Z, Li Z, You Q, Lu N, Guo Q. VI-14, a novel flavonoid derivative, inhibits migration and invasion of human breast cancer cells. *Toxicol Appl Pharmacol*. 2012 Jun 1;261(2):217-26.
- Chen FH, Lu N, Zhang HW, Zhao L, He LC, Sun HP, You QD, Li ZY, Guo QL. LYG-202 augments tumor necrosis factor- α -induced apoptosis via attenuating casein kinase 2-dependent nuclear factor- κ B pathway in HepG2 cells. *Mol Pharmacol*. 2012 Nov;82(5):958-71.
- Yao J, Hu R, Sun J, Lin B, Zhao L, Sha Y, Zhu B, You QD, Yan T, Guo QL. Oroxlylin A prevents inflammation-related tumor through down-regulation of inflammatory gene expression by inhibiting NF- κ B signaling. *Mol Carcinog*. 2014 Feb;53(2):145-58.
- Zhao J, Liu K, Lu J, Ma J, Zhang X, Jiang Y, Yang H, Jin G, Zhao G, Zhao M, Dong Z. Alternariol induces DNA polymerase β expression through the PKA-CREB signaling pathway. *Int J Oncol*. 2012 Jun;40(6):1923-8.
- Song X, Yao J, Wang F, Zhou M, Zhou Y, Wang H, Wei L, Zhao L, Li Z, Lu N, Guo Q. Wogonin inhibits tumor angiogenesis via degradation of

- HIF-1 α protein. *Toxicol Appl Pharmacol*. 2013 Sep 1;271(2):144-55.
16. Zhu L, Zhao L, Wang H, Wang Y, Pan D, Yao J, Li Z, Wu G, Guo Q. Oroxylin A reverses P-glycoprotein-mediated multidrug resistance of MCF7/ADR cells by G2/M arrest. *Toxicol Lett*. 2013 May 23; 219(2):107-15.
 17. Wu WY, Wang WY, Ma YL, Yan H, Wang XB, Qin YL, Su M, Chen T, Wang YP. Sodium tanshinone IIA silicate inhibits oxygen-glucose deprivation/recovery-induced cardiomyocyte apoptosis via suppression of the NF- κ B/TNF- α pathway. *Br J Pharmacol*. 2013 Jul;169(5):1058-71.
 18. Sun J, Fan L, Li M, Zhang Y, Cheng N. Decreasing Pin1 suppresses telomerase activity by NF- κ B in HCT116 cells colorectal carcinoma. *The Chinese-German Journal of Clinical Oncology*. 2013 April;12(4):181-187.
 19. Li K, Dan Z, Hu X, Ouzhu M, Ciren Y, Wang Z, Wang J, Yang X, Ze Y. CD14 overexpression upregulates TNF- α -mediated inflammatory responses and suppresses the malignancy of gastric carcinoma cells. *Mol Cell Biochem*. 2013 Apr;376(1-2):137-43.
 20. Yu Q, Shen XH, Li Y, Li RJ, Li J, Luo YY, Liu SF, Deng MY, Pei MF, Zhang GS. An intron mutation in the ACVRL1 may be associated with a transcriptional regulation defect in a Chinese family with hereditary hemorrhagic telangiectasia. *PLoS One*. 2013;8(2):e58031.
 21. Yan S, Fu Q, Zhou Y, Zhang N, Zhou Q, Wang X, Yuan Z, Wang X, Du J, Zhang J, Zhan L. Establishment of stable reporter expression for in vivo imaging of nuclear factor- κ B activation in mouse liver. *Theranostics*. 2013 Oct 15;3(11):841-50.
 22. Fang L, Xu Z, Wang GS, Ji FY, Mei CX, Liu J, Wu GM. Directed evolution of an LBP/CD14 inhibitory peptide and its anti-endotoxin activity. *PLoS One*. 2014 Jul 15;9(7):e101406.
 23. Wang G, Xiu P, Li F, Xin C, Li K. Vitamin A supplementation alleviates extrahepatic cholestasis liver injury through Nrf2 activation. *Oxid Med Cell Longev*. 2014;2014:273692. doi: 10.1155/2014/273692. Epub 2014 Jul 9.
 24. Zhao K, Song X, Huang Y, Yao J, Zhou M, Li Z, You Q, Guo Q, Lu N. Wogonin inhibits LPS-induced tumor angiogenesis via suppressing PI3K/Akt/NF- κ B signaling. *Eur J Pharmacol*. 2014 Aug 15;737:57-69.
 25. Yao LL, Liao CH, Huang G, Zhou Y, Rigali S, Zhang B, Ye BC. GlnR-mediated regulation of nitrogen metabolism in the actinomycete *Saccharopolyspora erythraea*. *Appl Microbiol Biotechnol*. 2014 Sep; 98(18):7935-48.
 26. Liao C, Rigali S, Cassani CL, Marcellin E, Nielsen LK, Ye BC. Control of chitin and N-acetylglucosamine utilization in *Saccharopolyspora erythraea*. *Microbiology*. 2014 Sep;160(Pt 9):1914-28.
 27. Xu X, Zhang Y, Li W, Miao H, Zhang H, Zhou Y, Li Z, You Q, Guo Q, Zhao L. Wogonin Reverses Multi-Drug Resistance of Human myelogenous leukemia K562/A02 cells via Downregulation of MRP1 Expression by Inhibiting Nrf2/ARE Signaling Pathway. *Biochem Pharmacol*. 2014 Nov 15;92(2):220-34.
 28. Li X, Jiang J, Shi S, Bligh SW, Li Y, Jiang Y, Huang D, Ke Y, Wang S. A RG-II type polysaccharide purified from *Aconitum coreanum* alleviates lipopolysaccharide-induced inflammation by inhibiting the NF- κ B signal pathway. *PLoS One*. 2014 Jun 13;9(6):e99697.
 29. Li X, Miao H, Zhang Y, Li W, Li Z, Zhou Y, Zhao L, Guo Q. Bone marrow microenvironment confers imatinib resistance to chronic myelogenous leukemia and oxorxylin A reverses the resistance by suppressing Stat3 pathway. *Arch Toxicol*. 2015 Jan;89(1):121-36.
 30. Fang Y, Chai Z, Wang D, Kuang T, Wu W, Lou W. DNA-PKcs deficiency sensitizes the human hepatoma HepG2 cells to cisplatin and 5-fluorouracil through suppression of the PI3K/Akt/NF- κ B pathway. *Mol Cell Biochem*. 2015 Jan;399(1-2):269-78.
 31. Zhao L, Miao HC, Li WJ, Sun Y, Huang SL, Li ZY, Guo QL. LW-213 induces G2/M cell cycle arrest through AKT/GSK3 β -catenin signaling pathway in human breast cancer cells. *Mol Carcinog*. 2016 May;55(5):778-92.
 32. Zhao Y, Wang X, Sun Y, Zhou Y, Yin Y, Ding Y, Li Z, Guo Q, Lu N. LYG-202 exerts antitumor effect on PI3K/Akt signaling pathway in human breast cancer cells. *Apoptosis*. 2015 Sep;20(9):1253-69.
 33. Wang W, Liu X, Wang X, Dong H, Ma C, Wang J, Liu B, Mao Y, Wang Y, Li T, Yang C, Guo Y. Structural and Functional Diversity of Nairovirus-Encoded Nucleoproteins. *J Virol*. 2015 Dec 1;89(23):11740-9.
 34. Dai L, Xu Y, Yu W, Liu S, Gao Y, Zhang L, Yuan B, Chen J, Ma T, Zhang J. Naturally occurring genetic mutations in the 5'-upstream regulatory region of bovine FSHB generate a novel cis-regulatory element that affects its expression. *Anim Genet*. 2015 Dec;46(6):693-6.
 35. Yang Y, Wu HJ, Lin L, Zhu QQ, Borriss R, Gao XW. A plasmid-born Rap-Phr system regulates surfactin production, sporulation and genetic competence in the heterologous host, *Bacillus subtilis* OKB105. *Appl Microbiol Biotechnol*. 2015 Sep;99(17):7241-5.
 36. Zhao L, Li W, Zhou Y, Zhang Y, Huang S, Xu X, Li Z, Guo Q. The overexpression and nuclear translocation of Trx-1 during hypoxia confers on HepG2 cells resistance to DDP, and GL-V9 reverses the resistance by suppressing the Trx-1/Ref-1 axis. *Free Radic Biol Med*. 2015 May;82:29-41.
 37. Liao CH, Yao L, Xu Y, Liu WB, Zhou Y, Ye BC. Nitrogen regulator GlnR controls uptake and utilization of non-phosphotransferase-system carbon sources in actinomycetes. *Proc Natl Acad Sci U S A*. 2015 Dec 22;112(51):15630-5.
 38. Zhang Y, Gao T, Hu S, Lin B, Yan D, Xu Z, Zhang Z, Mao Y, Mao H, Wang L, Wang G, Xiong Y, Zuo B. The Functional SNPs in the 5' Regulatory Region of the Porcine PPAR α Gene Have Significant Association with Fat Deposition Traits. *PLoS One*. 2015 Nov 24;10(11):e0143734.
 39. Yang L, Ni L, Duan Q, Wang X, Chen C, Chen S, Chaugai S, Zeldin DC, Tang JR, Wang DW. CYP epoxygenase 2J2 prevents cardiac fibrosis by suppression of transmission of pro-inflammation from cardiomyocytes to macrophages. *Prostaglandins Other Lipid Mediat*. 2015 Jan-Mar;116-117:64-75.
 40. Guo Y, Wang W, Sun Y, Ma C, Wang X, Wang X, Liu P, Shen S, Li B, Lin J, Deng F, Wang H, Lou Z. Crystal structure of the core region of hantavirus nucleocapsid protein reveals the mechanism for ribonucleoprotein complex formation. *J Virol*. 2015 Nov 11;90(2):1048-61.
 41. Zhang Y, Li W, Zhu M, Li Y, Xu Z, Zuo B. FHL3 differentially regulates the expression of MyHC isoforms through interactions with MyoD and pCREB. *Cell Signal*. 2016 Jan;28(1):60-73.
 42. Xia X, Yan C, Wu W, Zhou Y, Hou L, Zuo B, Xu D, Ren Z, Xiong Y. Characterization of the porcine peptidylarginine deiminase type VI gene (PADI6) promoter: Sp1 regulates basal transcription of the porcine PADI6. *Gene*. 2016 Jan 10;575(2 Pt 2):551-8.
 43. Wei L, Zhou Y, Qiao C, Ni T, Li Z, You Q, Guo Q, Lu N. Oroxylin A inhibits glycolysis-dependent proliferation of human breast cancer via promoting SIRT3-mediated SOD2 transcription and HIF1 α destabilization. *Cell Death Dis*. 2015 Apr 9;6:e1714.
 44. Liao CH, Xu Y, Rigali S, Ye BC. DasR is a pleiotropic regulator required for antibiotic production, pigment biosynthesis, and morphological development in *Saccharopolyspora erythraea*. *Appl Microbiol Biotechnol*. 2015 Dec;99(23):10215-24.
 45. Yao LL, Ye BC. The reciprocal regulation of GlnR and PhoP in response to nitrogen and phosphate limitations in *Saccharopolyspora erythraea*. *Appl Environ Microbiol*. 2015 Oct 30;82(1):409-20.
 46. Wang S, Liu K, Xiao L, Yang L, Li H, Zhang F, Lei L, Li S, Feng X, Li A, He J. Characterization of a novel DNA glycosylase from *S. sahachiroi* involved in the reduction and repair of azinomycin B induced DNA damage. *Nucleic Acids Res*. 2016 Jan 8;44(1):187-97.
 47. Jiang Q, Liu Z, Zhou Z, Wang L, Wang L, Yue F, Wang J, Wang H, Song L. Transcriptional activation and translocation of ancient NOS during immune response. *FASEB J*. 2016 Oct;30(10):3527-3540.
 48. Dong Y, Wang S, Chen J, Zhang Q, Liu Y, You C, Monroig Ó, Tocher DR, Li Y. Hepatocyte Nuclear Factor 4 α (HNF4 α) Is a Transcription Factor of Vertebrate Fatty Acyl Desaturase Gene as Identified in Marine Teleost *Siganus canaliculatus*. *PLoS One*. 2016 Jul 29;11(7):e0160361.
 49. You M, Jin J, Liu Q, Xu Q, Shi J, Hou Y. PPAR α Promotes Cancer Cell Glut1 Transcription Repression. *J Cell Biochem*. 2016 Dec 5. doi: 10.1002/jcb.25817. [Epub ahead of print]
 50. Wu XB, Liu Y, Wang GH, Xu X, Cai Y, Wang HY, Li YQ, Meng HF, Dai F, Jin JD. Mesenchymal stem cells promote colorectal cancer progression through AMPK/mTOR-mediated NF- κ B activation. *Sci Rep*. 2016 Feb 19;6:21420.
 51. Zhang Y, Li W, Zhu M, Li Y, Xu Z, Zuo B. FHL3 differentially regulates the expression of MyHC isoforms through interactions with MyoD and pCREB. *Cell Signal*. 2016 Jan;28(1):60-73.
 52. Xu X, Zhang X, Liu Y, Yang L, Huang S, Lu L, Wang S, Guo Q, Zhao L. BM microenvironmental protection of CML cells from imatinib through Stat5/NF- κ B signaling and reversal by Wogonin. *Oncotarget*. 2016 Apr 26;7(17):24436-54.
 53. Hong Y, Zhang H, Huang L, Li D, Song F. Overexpression of a Stress-Responsive NAC Transcription Factor Gene ONAC022 Improves

- Drought and Salt Tolerance in Rice. *Front Plant Sci.* 2016 Jan 22;7:4.
54. Liu X, Qin Z, Li X, Ma X, Gao B, Zhang Z. NF1, Sp1 and HSF1 are synergistically involved in sulfide-induced *sqr* activation in echiuran worm *Urechis unicinctus*. *Aquat Toxicol.* 2016 Jun;175:232-40.
55. Wu L, Guo H, Sun H, Zhang W, Sun C, Wang J. UNC119 mediates gambogic acid-induced cell-cycle dysregulation through the Gsk3 β / β -catenin pathway in hepatocellular carcinoma cells. *Anticancer Drugs.* 2016 Nov;27(10):988-1000.
56. Xia X, Yan C, Wu W, Zhou Y, Hou L, Zuo B, Xu D, Ren Z, Xiong Y. Characterization of the porcine peptidylarginine deiminase type VI gene (PADI6) promoter: Sp1 regulates basal transcription of the porcine PADI6. *Gene.* 2016 Jan 10;575(2 Pt 2):551-8.
57. Xu Y, Liao CH, Yao LL, Ye X, Ye BC. GlnR and PhoP directly regulate the transcription of genes encoding starch-degrading, amylolytic enzymes in *Saccharopolyspora erythraea*. *Appl Environ Microbiol.* 2016 Sep 16. pii: AEM.02117-16
58. Liu X, Zhang Z, Ma X, Li X, Zhou D, Gao B, Bai Y. Sulfide exposure results in enhanced *sqr* transcription through upregulating the expression and activation of HSF1 in echiuran worm *Urechis unicinctus*. *Aquat Toxicol.* 2016 Jan;170:229-39.
59. Zhuang K, Zhang J, Xiong M, Wang X, Luo X, Han L, Meng Y, Zhang Y, Liao W, Liu S. CDK5 functions as a tumor promoter in human colorectal cancer via modulating the ERK5-AP-1 axis. *Cell Death Dis.* 2016 Oct 13;7(10):e2415.
60. Wang S, Liu K, Xiao L, Yang L, Li H, Zhang F, Lei L, Li S, Feng X, Li A, He J. Characterization of a novel DNA glycosylase from *S. sahachiroi* involved in the reduction and repair of azinomycin B induced DNA damage. *Nucleic Acids Res.* 2016 Jan 8;44(1):187-97.

Version 2017.10.20