碧云天生物技术/Beyotime Biotechnology
订货热线：400-168-3301或800-8283301
订货e-mail：order@beyotime.com
技术咨询：info@beyotime.com
网址：http://www.beyotime.com

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- Zhao Y, Liu H, Li Y, Wu J, Greenlee AR, Yang C, Jiang Y. The role of miR-506 in transformed 16HBE cells induced by anti-benzo[a]pyrene-trans-7,8-dihydrodiol-9,10-epoxide. *Toxicol Lett.* 2011 Sep 10; 205(3):320-6.
- Cao Y, Mao X, Sun C, Zheng P, Gao J, Wang X, Min D, Sun H, Xie N, Cai J. Baicalin attenuates global cerebral ischemia/reperfusion injury in gerbils via anti-oxidative and anti-apoptotic pathways. *Brain Res Bull.* 2011 Jul 15;85(6):396-402.
- Duan J, Zhang Q, Zhao H, Du J, Bai F, Bai G. Cloning, expression, characterization and application of atcA, atcB and atcC from *Pseudomonassp.* for the production of L-cysteine. *Biotechnol Lett.* 2012 Jun;34(6):1101-6.
- Zhou Z , Chen S, Zhao H, Wang C, Gao K, Guo Y, Shen Z, Wang Y, Wang H, Mei X. Producol inhibits neural cell apoptosis via inhibition of mTOR signaling pathway after spinal cord injury. *Neuroscience.* 2016 Aug 4;329:193-200.
- Huang D, Li Y, Cui F, Chen J, Sun J. Purification and characterization of a novel polysaccharide-peptide complex from *Clinacanthus nutans*Lindau leaves. *Carbohydr Polym.* 2016 Feb 10;137:701-8.
- Hu J, Wang G, Zhao W, Gao W. In situ growth of a C-terminal interferon-alpha conjugate of a phospholipid polymer that outperforms PEGASYS in cancer therapy. *J Control Release.* 2016 Sep 10;237:71-7.
- Qiu Q, Jiang J, Lin L, Cheng S, Xin D, Jiang W, Shen J, Hu Z. Downregulation of RSK2 influences the biological activities of human osteosarcoma cells throughactivating AKT/mTOR signaling pathways. *Int J Oncol.* 2016 Jun;48(6):2508-20.
- Liu X, Xiao XY, Shou QY, Yan JF, Chen L, Fu HY, Wang JC. Bufalin inhibits pancreatic cancer by inducing cell cycle arrest via the c-Myc/NF-κB pathway. *J Ethnopharmacol.* 2016 Dec 4;193:538-545.
- Zhang M, Du Y, Lu R, Shu Y, Zhao W, Li Z, Zhang Y, Liu R, Yang T, Luo S, Gao M, Zhang Y, Zhang G, Liu J, Lu Y. Cholesterol Retards Senescence in Bone Marrow Mesenchymal Stem Cells By Modulating Autophagy andROS/p53/p21Cip1/Waf1 Pathway. *Oxid Med Cell Longev.* 2016;2016:7524308.
- Gao Y, Yang H, Fan Y, Li L, Fang J, Yang W. Hydrogen-Rich Saline Attenuates Cardiac and Hepatic Injury in Doxorubicin Rat Model by Inhibiting Inflammation and Apoptosis. *Mediators Inflamm.* 2016; 2016:1320365.
- Ouyang F, Huang H, Zhang M, Chen M, Huang H, Huang F, Zhou S. HMGB1 induces apoptosis and EMT in association with increased autophagy following H/R injury in cardiomyocytes. *Int J Mol Med.* 2016 Mar;37(3):679-89.
- Qin Z, Zhao L, Hu H, Jiang H, Yu L, Zeng S. Utilizing single- and double-transfected cell models expressing human organic anion transporter 1 and human cytochrome P450 1A2 to investigate the interactions with ingredients of herbal medicines. *Xenobiotica.* 2016 Aug

13. Liu L, Wang C, Lin Y, Xi Y, Li H, Shi S, Li H, Zhang W, Zhao Y, Tian Y, Xu C, Wang L. Suppression of calcium sensing receptor ameliorates cardiac hypertrophy through inhibition of autophagy. Mol Med Rep. 2016 Jul;14(1):111-20.
14. Wang M, Lu X. Exploring the Synergy between Cellobiose Dehydrogenase from Phanerochaete chrysosporium and Cellulase from Trichoderma reesei. Front Microbiol. 2016 Apr 29;7:620.
15. Wang C, Li Z, Han H, Luo G, Zhou B, Wang S, Wang J. Impairment of object recognition memory by maternal bisphenol A exposure is associated with inhibition of Akt and ERK/CREB/BDNF pathway in the male offspring hippocampus. Toxicology. 2016 Feb 3;341-343:56-64.
16. Chen Y, Hu Y, Liu S, Zheng H, Wu X, Huang Z, Li H, Peng B, Long J, Pan B, Huang C, Dong Q. Whole-body aerosol exposure of cadmium chloride (CdCl₂) and tetrabromobisphenol A (TBBPA) induced hepatic changes in CD-1 male mice. J Hazard Mater. 2016 Nov 15;318:109-16.
17. Wang X, Liu B, Dou Y, Fan H, Wang S, Li T, Ding C, Yu S. The Riemerella anatipestifer AS87_01735 Gene Encodes Nicotinamidase PncA, an Important VirulenceFactor. Appl Environ Microbiol. 2016 Sep 16;82(19):5815-23.
18. Liao ZJ, Liang RS, Shi SS, Wang CH, Yang WZ. Effect of baicalin on hippocampal damage in kainic acid-induced epileptic mice. Exp Ther Med. 2016 Sep;12(3):1405-1411.
19. Gao S, Yu HN, Wu YF, Liu XY, Cheng AX, Lou HX. Cloning and functional characterization of a phenolic acid decarboxylase from the liverwort *Conocephalumjaponicum*. Biochem Biophys Res Commun. 2016 Dec 9;481(3-4):239-244.
20. Li Y, Liu W, Li L, Hölscher C. Neuroprotective effects of a GIP analogue in the MPTP Parkinson's disease mouse model. Neuropharmacology. 2016 Feb;101:255-63.
21. Zhang D, Liu H, Zeng J, Miao X, Huang W, Chen H, Huang Y, Li Y, Ye D. Glucocorticoid exposure in early placentation induces preeclampsia in rats via interfering trophoblast development. Gen Comp Endocrinol. 2016 Jan 1;225:61-70.
22. Meng H, Song Y, Zhu J, Liu Q, Lu P, Ye N, Zhang Z, Pang Y, Qi J, Wu H. LRG1 promotes angiogenesis through upregulating the TGF β 1 pathway in ischemic rat brain. Mol Med Rep. 2016 Dec; 14(6):5535-5543.
23. Xiang S, Zeng Y, Xiong B, Qin Y, Huang X, Jiang Y, Luo W, Sooranna SR, Pinhu L. Transforming growth factor beta 1 induced endothelin-1 release is peroxisome proliferator-activatedreceptor gamma dependent in A549 cells. J Inflamm (Lond). 2016 Jun 10;13:19.

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