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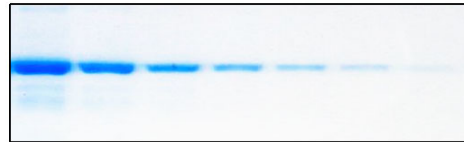
1. Liang W, Cheng J, Liu R, Wang JP, Mu JG, Wang QH, Wang HJ, Ma D. Peptide corresponding to the C terminus of tissue factor pathway inhibitor inhibits mesangial cell proliferation and activation in vivo. *Peptides*. 2009;30(12):2330-6.
2. Yan L, Xu G, Qiao T, Chen W, Yuan S, Li X. CpG-ODN 7909 Increases Radiation Sensitivity of Radiation-Resistant Human Lung Adenocarcinoma Cell Line by Overexpression of Toll-Like Receptor 9. *Cancer Biother Radiopharm*. 2013 Sep;28(7):559-64.
3. Zhang HH, Pang M, Dong W, Xin JX, Li YJ, Zhang ZC, Yu L, Wang PY, Li BS, Xie SY. miR-511 induces the apoptosis of radioresistant lung adenocarcinoma cells by triggering BAX. *Oncol Rep*. 2014 Mar;31(3):1473-9.
4. Wang X, Zhao B, Li X. Dexmedetomidine attenuates isoflurane-induced cognitive impairment through antioxidant, anti-inflammatory and anti-apoptosis in aging rat. *Int J Clin Exp Med*. 2015 Oct 15;8(10):17281-8.
5. Xing N, Qiao T, Zhuang X, Yuan S, Zhang Q, Xu G. CpG oligodeoxynucleotide 7909 enhances radiosensitivity via downregulating Oct-4 expression in radioresistant lung cancer cells. *Onco Targets Ther*. 2015 Jun 12;8:1443-9.
6. Zhuang X, Qiao T, Xu G, Yuan S, Zhang Q, Chen X. Combination of nadroparin with radiotherapy results in powerful synergistic antitumor effects in lung adenocarcinoma A549 cells. *Oncol Rep*. 2016 Oct;36(4):2200-6.
7. Ma LM, Liang ZR, Zhou KR, Zhou H, Qu LH. 27-Hydroxycholesterol increases Myc protein stability via suppressing PP2A, SCP1 and FBW7 transcription in MCF-7 breast cancer cells. *Biochem Biophys Res Commun*. 2016 Nov 18;480(3):328-333.
8. Zhang K, Zhu X, Lu Y. Effect of Mild Heating on Human Lens Epithelial Cells: A Possible Model of Lens Aging. *Sci Rep*. 2016 Oct 11;6:33917.

Version 2016.12.05