Anti-Flag Affinity Gel (Anti-Flag亲和凝胶)

产品编号	产品名称	包装
P2271-0.5ml	Anti-Flag Affinity Gel (Anti-Flag亲和凝胶)	0.5ml
P2271-2ml	Anti-Flag Affinity Gel (Anti-Flag亲和凝胶)	2ml
P2271-10ml	Anti-Flag Affinity Gel (Anti-Flag亲和凝胶)	10ml

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Characteristics	Description
Product content	50% settled gel in TBS with preservative
Matrix	4% cross - linked agarose
Average bead size	90 μ m
Antibody	Mouse monoclonal antibody against Flag-tag, clone G1
Isotype	IgG2b
M.W. of antibody	Approximately 150kDa
Antibody concentration	Approximately 8mg Flag antibody per ml settled gel
Binding capacity	Approximately 1mg Flag - tagged protein per ml settled gel
Gel reuse	The gel can be recycled for at least 4 times. If maintained properly, the gel can be reused 10 times with minimum loss of binding capability.
Elution method	Acid, alkaline, neutral, peptide competitive or SDS - PAGE loading buffer elutions. Note: If elute with SDS-PAGE loading buffer, the light (~25kDa) and heavy (~50kDa) chain of G1 antibody will be denatured and release from the gel.
Reagents compatibility	Compatible with commonly used reagents at certain concentrations
Application	Protein purification、Pull down、IP、Co-IP

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Reagent	Maximum tolerable concentration	Comments
EDTA	5mM	Higher concentration of chelating agent will reduce purification efficiency with less target protein recovery.
β - ME、 DTT	100mM for β - ME; 80mM for DTT	Reducing agents will reduce disulfide bonds in the G1 antibody on the gel. Avoid reducing agents or keep at low concentration (<10mM) during purification process. When the reducing agents reach the maximum tolerable concentration mentioned here, there will be bands of antibody heavy chain (~50kDa) and light chain (~25kDa) in SDS - PAGE analysis. The gel CANNOT BE REUSED if samples containing higher concentration of reducing reagents are applied to the gel.
Tween 20、 Triton X - 100	5% for Tween 20 and Triton X - 100	The concentration of Tween 20 and Triton X - 100 should not exceed 5%.
SDS	Not suggested	SDS will denature the G1 antibody on the gel.
NP - 40	4%	Higher concentration of NP - 40 will reduce purification efficiency with less target protein recovery.
GuHCl、Urea	0.3M for GuHCl; 1.5M for Urea	Chaotropic reagents will denature the target Flag - tagged protein. Do not exceed 0.3M GuHCl or 1.5M Urea.
Glycerol	20%	Higher concentration of glycerol will interfere with the binding of Flag tagged protein.
NaCl	0.5M	Higher concentration of NaCl will reduce purification efficiency with less target protein recovery.

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Problem	Possible Causes	Solution
Large amount of Flag-tagged protein found in the flow through.	Binding time is not enough.	If using batch method, increase the binding time experimentally; If using column method, use a lower flow rate when loading samples.
	Column is overloaded.	Reduce the amount of the sample added to the gel or increase the amount of gel.
	Flag - tag is not accessible to gel.	Expose the epitope tag by adding low amount of denaturant to the protein extract (dialysis may be needed before applying onto gel), or fuse Flag tag to the other terminus of the target protein.
	Gel has not been regenerated since last purification.	Perform gel regeneration procedure prior to binding.
	Reagent compatibility problem.	Dialyze the sample against TBS before purification procedure.
	The target protein has been degraded.	1. Prepare fresh lysates. Avoid using frozen lysates. 2. Perform purification at lower temperature, such as 4°C. 3. Use appropriate protease inhibitors in the lysate or increase their concentrations to prevent degradation of the Flag fusion protein.
Very few or no Flag - tagged protein exists in the eluate.	Protein is not completely eluted.	Change elution methods.
	No target protein expressed.	Make sure the protein of interest contains the Flag-tag by Western blot or dot blot analyses.
	Very low protein expression level.	1. Use larger volume of cell lysate. 2. Optimize expression conditions to raise the protein expression level.
	Washes are too stringent.	1. Reduce the number of washes. 2. Avoid adding high concentrations of NaCl to the mixture. 3. Use solutions that contain less or no detergent
	Incubation times are inadequate.	Increase the incubation times with the affinity gel (from several hours to overnight).
	Interfering substance is present in sample.	1. Lysates containing high concentrations of DTT, 2-mercaptoethanol, or other reducing agents may destroy antibody function, and must be avoided. 2. Excessive detergent concentrations may interfere with the antibody-antigen interaction. Detergent levels in buffers may be reduced by dilution.
	Detection system is inadequate.	If Western blot detection is used: 1. Check primary and secondary antibodies using proper controls to confirm binding and reactivity. 2. Verify that the transfer was adequate by staining the membrane with Ponceau S. 3. Use fresh detection substrate or try a different detection system.
Multiple protein bands found in the eluate.	The protein is not stable at room temperature.	Purify the target protein at lower temperature, such as 4°C.
	Protein degradation due to proteases activity during purification process.	Add protease inhibitors to cell lysate.
	Non - specific binding.	1. Prepare cell lysate again. 2. Add additional wash steps.
Background is too high.	Proteins bind nonspecifically to the Anti-Flag monoclonal antibody, the gel beads, or the microcentrifuge tubes.	1. Pre-clear lysate with Mouse IgG-Agarose (P2055+A7028) to remove nonspecific binding proteins. 2. After suspending beads for the final wash, transfer entire sample to a clean microcentrifuge tube before centrifugation.

	Washes are insufficient.	<ol style="list-style-type: none"> 1. Increase the number of washes. 2. Prolong duration of the washes, incubating each wash for at least 15 minutes. 3. Increase the salt and/or detergent concentrations in the wash solutions. 4. Centrifuge at lower speed to avoid nonspecific trapping of denatured proteins from the lysate during the initial centrifugation of the affinity gel complexes.
--	--------------------------	---

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产品编号	产品名称	包装
P2271-0.5ml	Anti-Flag Affinity Gel (Anti-Flag亲和凝胶)	0.5ml
P2271-2ml	Anti-Flag Affinity Gel (Anti-Flag亲和凝胶)	2ml
P2271-10ml	Anti-Flag Affinity Gel (Anti-Flag亲和凝胶)	10ml
P9801-1mg	3X Flag Peptide (3X Flag多肽)	5mg/ml×0.2ml
P9801-5mg	3X Flag Peptide (3X Flag多肽)	5mg/ml×1ml
P9801-25mg	3X Flag Peptide (3X Flag多肽)	25mg

Version 2021.05.13