碧云天生物技术/Beyotime Biotechnology
订货热线: 400-168-3301或800-8283301
订货e-mail: order@beyotime.com
技术咨询: info@beyotime.com
网址: http://www.beyotime.com

SNP (NO供体)

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1. Wu W, Xu X, Dai Y, Xia L. Therapeutic effect of the saponin fraction from Clematis chinensis Osbeck roots on osteoarthritis induced by monosodium iodoacetate through protecting articular cartilage. *Phytother Res.* 2010 Apr;24(4):538-46.
2. Fu YB, Yin HY, Wang WX, Wang MY, Zhang HY, Zhao XM, Du YG. β -1,3-Glucan with different degree of polymerization induced different defense responses in tobacco Carbohydrate Polymers. 2011 Aug 15;86(2):774-782.
3. Zhang Y, He LS, Zhang G, Xu Y, Lee OO, Matsumura K, Qian PY. The regulatory role of the NO/cGMP signal transduction cascade during larval attachment and metamorphosis of the barnacle Balanus (=Amphibalanus) amphitrite. *J Exp Biol.* 2012 Nov 1;215(Pt 21):3813-22.
4. He Z, Tan Z, Sun Z, Beauchemin KA, Tang S, Zhou C, Han X, Wang M, Wu D. Unchanged interleukin 6 level of protein and energy restricted goats during late gestation: the role of elevated blood nitric oxide. *J Endocrinol.* 2012 Apr;213(1):59-65.
5. Huang X, Zhou Y, Ma J, Wang N, Zhang Z, Ji J, Ding Q, Chen G. Nitric oxide mediated effects on reproductive toxicity caused by carbon disulfide in male rats. *Environ Toxicol Pharmacol.* 2012 Nov;34(3):679-87.
6. Wu W, Gao X, Xu X, Luo Y, Liu M, Xia Y, Dai Y. Saponin-rich fraction from Clematis chinensis Osbeck roots protects rabbit chondrocytes against nitric oxide-induced apoptosis via preventing mitochondria impairment and caspase-3 activation. *Cytotechnology.* 2013 Mar;65(2):287-95.
7. Sun GG, Hu WN, Zhang J, Li CL, Yang CR. Effect of nitric oxide on esophageal cancer cell line TE-1. *Chin Med Sci J.* 2013 Mar;28(1):44-9.
8. Xiao L, Liu W, Li J, Xie Y, He M, Fu J, Jin W, Shao C. Irradiated U937 cells trigger inflammatory bystander responses in human umbilical vein endothelial cells through the p38 pathway. *Radiat Res.* 2014 Jul;182(1):111-21.
9. Zhang G, Wong YH, Zhang Y, He LS, Xu Y, Qian PY. Nitric oxide inhibits larval settlement in Amphibalanus amphitrite cyprids by repressing muscle locomotion and molting. *Proteomics.* 2015 Nov;15(22):3854-64.
10. Zhou L, Zhang H, Wu J. Effects of nitric oxide on the biological behavior of HepG2 human hepatocellular carcinoma cells. *Exp Ther Med.* 2016 May;11(5):1875-1880.
11. Shen J, Tian Y, Li Y, Ma R, Zhang Q, Zhang J, Fang J. Bactericidal Effects against S. aureus and Physicochemical Properties of Plasma Activated Water stored at different temperatures. *Sci Rep.* 2016 Jun 27;6:28505.
12. Wu KW, Kou ZW, Mo JL, Deng XX, Sun FY. Neurovascular coupling protects neurons against hypoxic injury via inhibition of potassium currents by generation of nitric oxide in direct neuron and endothelium cocultures. *Neuroscience.* 2016 Oct 15;334:275-282.
13. Ji YY, Ma YJ, Wang JW. Cytoprotective role of nitric oxide in HepG2 cell apoptosis induced by hypocrellin B photodynamic treatment. *J Photochem Photobiol B.* 2016 Oct;163:366-73.
14. Zhai L, Xiao D, Sun C, Wu T, Han Z, Zhang X, Xu X, Wang Y. Nitric oxide signaling is involved in the response to iron deficiency in the woody plant Malus xiaojinensis. *Plant Physiol Biochem.* 2016 Dec;109:515-524.

15. Zhang G, Yan GY, Yang XX, Wong YH, Sun J, Zhang Y, He LS, Xu Y, Qian PY. Characterization of Arginine Kinase in the Barnacle Amphibalanus Amphitrite and Its Role in the Larval Settlement. *J Exp Zool B Mol Dev Evol.* 2016 Jun;326(4):237-49.
16. Zhao W, Wu C, Chen X. Cryptotanshinone inhibits oxidized LDL-induced adhesion molecule expression via ROS dependent NF- κ B pathways. *Cell Adh Migr.* 2016 May 3;10(3):248-58.
17. Guo Y, Deng Y, Huang Z, Luo Q, Peng Y, Chen J, Jiang H, Ye J, Li J. EBP50 induces apoptosis in macrophages by upregulating nitric oxide production to eliminate intracellular Mycobacterium tuberculosis. *Sci Rep.* 2016 Jan 5;6:18961.

Version 2016.12.08