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1. Wang DW, Hu YL, Liu YJ, Lin ZP. Studies on the Function of Vitreoscilla Hemoglobin Gene in Improving Sur factin Producing Rate of Bacillus subtilis. Biotechnology bulletin. 2007;5:116-9.
2. He Z, Sun X, Mei G, Yu S, Li N. Nonclassical secretion of human catalase on the surface of CHO cells is more efficient than classical secretion. Cell Biol Int. 2008 Apr;32(4):367-73.
3. He Z, Yu S, Mei G, Zheng M, Wang M, Dai Y, Tang B, Li N. Maternally transmitted milk containing recombinant human catalase provides protection against oxidation for mouse offspring during lactation. Free Radic Biol Med. 2008 Oct 15;45(8):1135-42.
4. Zhu M, Gu F, Shi J, Hu J, Hu Y, Zhao Z. Increased oxidative stress and astrogliosis responses in conditional double-knockout mice of Alzheimer-like presenilin-1 and presenilin-2. Free Radic Biol Med. 2008 Nov 15;45(10):1493-9.
5. Wang XZ, Liu SS, Sun Y, Wu JY, Zhou YL, Zhang JH. Beta-cypermethrin impairs reproductive function in male mice by inducing oxidative stress. Theriogenology. 2009 Sep 15;72(5):599-611.
6. Qian J, Jiang F, Wang B, Yu Y, Zhang X, Yin Z, Liu C. Ophiopogonin D prevents H₂O₂-induced injury in primary human umbilical vein endothelial cells. J Ethnopharmacol. 2010;128(2):438-45.
7. Huang X, Zhuang J, Teng X, Li L, Chen D, Yan X, Tang F. The promotion of human malignant melanoma growth by mesoporous silica nanoparticles through decreased reactive oxygen species. Biomaterials. 2010;31(24):6142-53.
8. Chen Y, Chang L, Li W, Rong Z, Liu W, Shan R, Pan R. Thioredoxin protects fetal type II epithelial cells from hyperoxia-induced injury. Pediatr Pulmonol. 2010 Dec;45(12):1192-200.
9. Liu XM, Zhou Q, Xu SZ, Wakamatsu K, Lei TC. Maintenance of immune hyporesponsiveness to melanosomal proteins by DHICA-mediated antioxidation: Possible implications for autoimmune vitiligo. Free Radic Biol Med. 2011 May 1;50(9):1177-85.
10. Zhou J, Li Y, Yan G, Bu Q, Lv L, Yang Y, Zhao J, Shao X, Deng Y, Zhu R, Zhao Y, Cen X. Protective role of taurine against morphine-induced neurotoxicity in C6 cells via inhibition of oxidative stress. Neurotox Res. 2011 Nov;20(4):334-42.
11. Min L, He S, Chen Q, Peng F, Peng H, Xie M. Comparative proteomic analysis of cellular response of human airway epithelial cells (A549) to benzo(a)pyrene. Toxicol Mech Methods. 2011 Jun;21(5):374-82.
12. Yang L, Zheng XL, Sun H, Zhong YJ, Wang Q, He HN, Shi XW, Zhou B, Li JK, Lin Y, Zhang L, Wang X. Catalase suppression-mediated H₂O₂ accumulation in cancer cells by wogonin effectively blocks tumor necrosis factor-induced NF-κB activation and sensitizes apoptosis.

- Cancer Sci. 2011 Apr;102(4):870-6.
13. Fan J, Yin JJ, Ning B, Wu X, Hu Y, Ferrari M, Anderson GJ, Wei J, Zhao Y, Nie G. Direct evidence for catalase and peroxidase activities of ferritin-platinum nanoparticles. *Biomaterials*. 2011 Feb;32(6):1611-8.
 14. Zhang Y, Fu D, Yu F, Liu Q, Yu Z. Two catalase homologs are involved in host protection against bacterial infection and oxidative stress in *Crassostrea hongkongensis*. *Fish Shellfish Immunol*. 2011 Dec;31(6):894-903.
 15. Li W, Wu Y, Ren C, Lu Y, Gao Y, Zheng X, Zhang C. The activity of recombinant human neuroglobin as an antioxidant and free radical scavenger. *Proteins*. 2011 Jan;79(1):115-25.
 16. Wang J, Sun P, Bao Y, Liu J, An L. Cytotoxicity of single-walled carbon nanotubes on PC12 cells. *Toxicol In Vitro*. 2011 Feb;25(1):242-50.
 17. Shi X, Gu A, Ji G, Li Y, Di J, Jin J, Hu F, Long Y, Xia Y, Lu C, Song L, Wang S, Wang X. Developmental toxicity of cypermethrin in embryo-larval stages of zebrafish. *Chemosphere*. 2011 Oct;85(6):1010-6.
 18. Wang S, Duan Y, Zeng Q, Zhao Z, Wang X. [D-Ala2, D-Leu5]-Enkephalin (Dadle) Reduces Inflammation Responses after Transient Global Ischemia *Advanced Materials Research*. 2011 Sep;345:343-48.
 19. Zhao G, Yao-Yue C, Qin GW, Guo LH. Luteolin from Purple Perilla mitigates ROS insult particularly in primary neurons. *Neurobiol Aging*. 2012 Jan;33(1):176-86.
 20. Deng X, Wu K, Wan J, Li L, Jiang R, Jia M, Jing Y, Zhang L. Aminotriazole attenuated carbon tetrachloride-induced oxidative liver injury in mice. *Food Chem Toxicol*. 2012 Jun 8;50(9):3073-8.
 21. Cao L, Li Z, Zhang R, Wu Y, Li W, Cao Z. StCT2, a new antibacterial peptide characterized from the venom of the scorpion *Scorpiopstibetanus*. *Peptides*. 2012 Aug;36(2):213-20.
 22. Wang H, Meng XL, Xu JP, Wang J, Wang H, Ma CW. Production, purification, and characterization of the cecropin from *Plutella xylostella*, pxCeCA1 using an intein-induced self-cleavable system in *Escherichia coli*. *Appl Microbiol Biotechnol*. 2012 May;94(4):1031-9.
 23. Shao X, Hu Z, Hu C, Bu Q, Yan G, Deng P, Lv L, Wu D, Deng Y, Zhao J, Zhu R, Li Y, Li H, Xu Y, Yang H, Zhao Y, Cen X. Taurine protects methamphetamine-induced developmental angiogenesis defect through antioxidant mechanism. *Toxicol Appl Pharmacol*. 2012 May 1;260(3):260-70.
 24. Li NS, Luo XJ, Dai Z, Liu B, Zhang YS, Yang ZC, Peng J. Beneficial effects of capsiate on ethanol-induced mucosal injury in rats are related to stimulation of calcitonin gene-related Peptide release. *Planta Med*. 2012 Jan;78(1):24-30.
 25. Liu Z, Li S, Cai Y, Wang A, He Q, Zheng C, Zhao T, Ding X, Zhou X. Manganese superoxide dismutase induces migration and invasion of tongue squamous cell carcinoma via H2O2-dependent Snail signaling. *Free Radic Biol Med*. 2012 Jul 1;53(1):44-50.
 26. Cao L, Dai C, Li Z, Fan Z, Song Y, Wu Y, Cao Z, Li W. Antibacterial activity and mechanism of a scorpion venom Peptide derivative in vitro and in vivo. *PLoS One*. 2012;7(7):e40135.
 27. Liu JH, Qi ZM, Huang Q, Wei XL, Ke ZG, Fang YS, Tian YC, Yu ZG. Study of energetic-particle-irradiation induced biological effect on *Rhizopus oryzae* through Synchrotron-FTIR micro-spectroscopy. *Journal of Molecular Structure*. 2012 Jan;1031(16):1-8.
 28. Li XC, Xing YZ, Jiang X, Qiao J, Tan HL, Zhou B. Identification And Characterization Of The Catalase Gene Pycat From The Red Alga *Pyropia Yezoensis* (Bangiales, Rhodophyta) 1. *Journal of Phycology*. 2012 Jun;48(3):664-9.
 29. Wang J, Sun P, Bao Y, Dou B, Song D, Li Y. Vitamin E renders protection to PC12 cells against oxidative damage and apoptosis induced by single-walled carbon nanotubes. *Toxicol In Vitro*. 2012 Feb;26(1):32-41.
 30. Sun WH, Liu F, Chen Y, Zhu YC. Hydrogen sulfide decreases the levels of ROS by inhibiting mitochondrial complex IV and increasing SOD activities in cardiomyocytes under ischemia/reperfusion. *Biochem Biophys Res Commun*. 2012 May 4;421(2):164-9.
 31. Zhou T, Zong R, Zhang Z, Zhu C, Pan F, Xiao X, Liu Z, He H, Ma JX, Liu Z, Zhou Y. SERPINA3K Protects against Oxidative Stress via Modulating ROS Generation/Degradation and KEAP1-NRF2 Pathway in the Corneal Epithelium. *Invest Ophthalmol Vis Sci*. 2012 Jul 27;53(8):5033-43.
 32. Tie L, Yang HQ, An Y, Liu SQ, Han J, Xu Y, Hu M, Li WD, Chen AF, Lin ZB, Li XJ. Ganoderma lucidum polysaccharide accelerates refractory wound healing by inhibition of mitochondrial oxidative stress in type 1 diabetes. *Cell Physiol Biochem*. 2012;29(3-4):583-94.
 33. Shi H, Wang Y, Cheng Z, Ye T, Chan Z. Analysis of natural variation in bermudagrass (*Cynodon dactylon*) reveals physiological responses underlying drought tolerance. *PLoS One*. 2012;7(12):e53422.
 34. Zheng J, Yang B, Yu Y, Chen Q, Huang T, Li D. Ganoderma lucidum polysaccharides exert anti-hyperglycemic effect on streptozotocin-induced diabetic rats through affecting β -cells. *Comb Chem High Throughput Screen*. 2012 Aug;15(7):542-50.
 35. Ren C, Li Y, Han R, Gao D, Li W, Shi J, Hoogewijs D, Braeckman BP, De Henau S, Lu Y, Qu W, Gao Y, Wu Y, Li Z, Liu H, Wang Z, Zhang C. GLB-13 is associated with oxidative stress resistance in *Caenorhabditis elegans*. *IUBMB Life*. 2013 May;65(5):423-34.
 36. Sun P, Chen JY, Li J, Sun MR, Mo WC, Liu KL, Meng YY, Liu Y, Wang F, He RQ, Hua Q. The protective effect of geniposide on human neuroblastoma cells in the presence of formaldehyde. *BMC Complement Altern Med*. 2013 Jul 1;13:152.
 37. Wang X, Zhang Y, Li T, Tian W, Zhang Q, Cheng Y. Generation 9 polyamidoamine dendrimer encapsulated platinum nanoparticle mimics catalase size, shape, and catalytic activity. *Langmuir*. 2013 Apr 30;29(17):5262-70.
 38. Wang M, Wang L, Zhou Z, Gao Y, Wang L, Shi X, Gai Y, Mu C, Song L. The molecular characterization of a catalase from Chinese mitten crab *Eriocheir sinensis*. *Int J Immunogenet*. 2013 Jun;40(3):230-40.
 39. Shi H, Ye T, Wang Y, Chan Z. Arabidopsis ALTERED MERISTEM PROGRAM 1 negatively modulates plant responses to abscisic acid and dehydration stress. *Plant Physiol Biochem*. 2013 Apr 3;67C:209-216.
 40. Zhang H, Shao D, Wu Y, Dai B, Cai C, Fang W, Ye B, Zhang Y, Liu J, Jia X. Regulation of nodularin-induced apoptosis by epigallocatechin-3-gallate on fish lymphocytes in vitro. *Fish Shellfish Immunol*. 2013 May;34(5):1085-93.
 41. Yu CH, Liu ZY, Sun W, Li YJ, Zhang DS, Pan RT, Sun ZL. Effect of Danofloxacin on Reactive Oxygen Species Production, Lipid Peroxidation and Antioxidant Enzyme Activities in Kidney Tubular Epithelial Cell Line, LLC-PK1. *Basic Clin Pharmacol Toxicol*. 2013 Dec;113(6):377-84.
 42. Chen Z, Wang D, Gu C, Liu X, Pei W, Li J, Cao Y, Jiao Y, Tong J, Nie J. Down-regulation of let-7 microRNA increased K-ras expression in lung damage induced by radon. *Environ Toxicol Pharmacol*. 2015 Sep;40(2):541-8.
 43. He Z, Zhang X, Chen C, Wen Z, Hoopes SL, Zeldin DC, Wang DW. Cardiomyocyte-specific expression of CYP2J2 prevents development of cardiac remodeling induced by angiotensin II. *Cardiovasc Res*. 2015 Mar 1;105(3):304-17.
 44. Luo C, Li Y, Yang L, Wang X, Long J, Liu J. Superparamagnetic iron oxide nanoparticles exacerbate the risks of reactive oxygen species-mediated external stresses. *Arch Toxicol*. 2015 Mar;89(3):357-69.
 45. Cai W, Liu W, Wang WS, Fu ZW, Han TT, Lu YT. Overexpression of Rat Neurons Nitric Oxide Synthase in Rice Enhances Drought and Salt Tolerance. *PLoS One*. 2015 Jun 29;10(6):e0131599.
 46. Hu Y, Li R, Yang H, Luo H, Chen Z. Sirtuin 6 is essential for sodium sulfide-mediated cytoprotective effect in ischemia/reperfusion-stimulated brain endothelial cells. *J Stroke Cerebrovasc Dis*. 2015 Mar;24(3):601-9.
 47. Chen L, Bi B, Zeng J, Zhou Y, Yang P, Guo Y, Zhu J, Yang Q, Zhu N, Liu T. Rosiglitazone ameliorates senescence-like phenotypes in a cellular photoaging model. *J Dermatol Sci*. 2015 Mar;77(3):173-81.
 48. Su H, Liu DD, Zhao M, Hu WL, Xue SS, Cao Q, Le XY, Ji LN, Mao ZW. Dual-Enzyme Characteristics of Polyvinylpyrrolidone-Capped Iridium Nanoparticles and Their Cellular Protective Effect against H2O2-Induced Oxidative Damage *ACS Appl Mater Interfaces*. 2015 Apr 22;7(15):8233-42.
 49. Keogh K, Waters SM, Kelly AK, Wylie AR, Kenny DA. Effect of feed restriction and subsequent re-alimentation on hormones and genes of the somatotrophic axis in cattle. *Physiol Genomics*. 2015 Jul;47(7):264-73.
 50. Chen W, Su H, Xu Y, Bao T, Zheng X. Protective effect of wild raspberry (*Rubus hirsutus* Thunb.) extract against acrylamide-induced oxidative damage is potentiated after simulated gastrointestinal digestion. *Food Chem*. 2016 Apr 1;196:943-52.
 51. Dou T, Yan M, Wang X, Lu W, Zhao L, Lou D, Wu C, Chang X, Zhou Z. Nrf2/ARE Pathway Involved in Oxidative Stress Induced by Paraquat in Human Neural Progenitor Cells. *Oxid Med Cell Longev*. 2016;2016:8923860.
 52. Sun H, Li G, An T, Zhao H, Wong PK. Unveiling the photoelectron catalytic inactivation mechanism of *Escherichia coli*: Convincing

- evidence from responses of parent and anti-oxidation single gene knockout mutants. *Water Res.* 2016 Jan 1;88:135-43.
53. Lin J, Zou Y, Cao K, Ma C, Chen Z. The impact of heterologous catalase expression and superoxide dismutase overexpression on enhancing the oxidative resistance in *Lactobacillus casei*. *J Ind Microbiol Biotechnol.* 2016 May;43(5):703-11.
 54. Dong Y, Cui P, Li Z, Zhang S. Aging asymmetry: systematic survey of changes in age-related biomarkers in the annual fish *Nothobranchius guentheri*. *Fish Physiol Biochem.* 2016 Sep 10.[Epub ahead of print]
 55. Meng J, Lv Z, Qiao X, Li X, Li Y, Zhang Y, Chen C. The decay of Redox-stress Response Capacity is a substantive characteristic of aging: Revising the redox theory of aging. *Redox Biol.* 2016 Dec 28;11:365-374.
 56. Huang H, Xiao X, Lin F, Grossart HP, Nie Z, Sun L, Xu C, Shi J. Continuous-release beads of natural allelochemicals for the long-term control of cyanobacterial growth: Preparation, release dynamics and inhibitory effects. *Water Res.* 2016 May 15;95:113-23.
 57. Wang J, Chen X, Wang F, Zhang J, Li P, Li Z, Xu J, Gao F, Jin C, Tian H, Zhang J, Li W, Lu L, Xu GT. OFD1, as a Ciliary Protein, Exhibits Neuroprotective Function in Photoreceptor Degeneration Models. *PLoS One.* 2016 May 19;11(5):e0155860.
 58. Yao C, Li G, Qian Y, Cai M, Yin H, Xiao L, Tang W, Guo F, Shi B. Protection of Pentoxifylline against Testis Injury Induced by Intermittent Hypobaric Hypoxia. *Oxid Med Cell Longev.* 2016;2016:3406802.
 59. Jiao Y, Ma S, Wang Y, Li J, Shan L, Liu Q, Liu Y, Song Q, Yu F, Yu H, Liu H, Huang L, Chen J. N-Acetyl Cysteine Depletes Reactive Oxygen Species and Prevents Dental Monomer-Induced Intrinsic Mitochondrial Apoptosis In Vitro in Human Dental Pulp Cells. *PLoS One.* 2016 Jan 25;11(1):e0147858.
 60. Wang M, Zhao X, Xiao Z, Yin X, Xing T, Xia G. A wheat superoxide dismutase gene TaSOD2 enhances salt resistance through modulating redox homeostasis by promoting NADPH oxidase activity. *Plant Mol Biol.* 2016 May;91(1-2):115-30.
 61. Ran B, Li M, Li Y, Lin Y, Liu W, Luo Q, Fu Y, Tang Q, Yang Y, Pu Y. Everolimus (RAD001) inhibits the proliferation of rat vascular smooth muscle cells by up-regulating the activity of the p27/kip1 gene promoter. *Anatol J Cardiol.* 2016 Jun;16(6):385-91.
 62. Han Y, Su J, Liu X, Zhao Y, Wang C, Li X. Naringin alleviates early brain injury after experimental subarachnoid hemorrhage by reducing oxidative stress and inhibiting apoptosis. *Brain Res Bull.* 2016 Dec 21. pii: S0361-9230(16)30479-8.
 63. Liu Z, Gan L, Chen Y, Luo D, Zhang Z, Cao W, Zhou Z, Lin X, Sun C. Mark4 promotes oxidative stress and inflammation via binding to PPAR γ and activating NF- κ B pathway in mice adipocytes. *Sci Rep.* 2016 Feb 18;6:21382.
 64. Chen W, Su H, Xu Y, Jin C. In vitro gastrointestinal digestion promotes the protective effect of blackberry extract against acrylamide-induced oxidative stress. *Sci Rep.* 2017 Jan 13;7:40514.
 65. Chen YQ, Liu XG, Zhao W, Cui H, Ruan J, Yuan Y, Tu Z. MET18 Deficiency Increases the Sensitivity of Yeast to Oxidative Stress and Shortens Replicative Lifespan by Inhibiting Catalase Activity. *Biomed Res Int.* 2017;2017:7587395.
 66. Dong Y, Cui P, Li Z, Zhang S. Aging asymmetry: systematic survey of changes in age-related biomarkers in the annual fish *Nothobranchius guentheri*. *Fish Physiol Biochem.* 2017 Apr;43(2):309-319.
 67. Chen W, Su H, Xu Y, Jin C. In vitro gastrointestinal digestion promotes the protective effect of blackberry extract against acrylamide-induced oxidative stress. *Sci Rep.* 2017 Jan 13;7:40514.
 68. Zhu Z, Fan X, Pan Y, Lu Y, Zeng W. Trehalose improves rabbit sperm quality during cryopreservation. *Cryobiology.* 2017 Apr;75:45-51.
 69. Zhang GM, Deng MT, Lei ZH, Wan YJ, Nie HT, Wang ZY, Fan YX, Wang F, Zhang YL. Effects of NRF1 on steroidogenesis and apoptosis in goat luteinized granulosa cells. *Reproduction.* 2017 Aug;154(2):111-122.
 70. Gao Y, Zhang J, Liu Y, Zhang S, Wang Y, Liu B, Liu H, Li R, Lv C, Song X. Regulation of TERRA on telomeric and mitochondrial functions in IPF pathogenesis. *BMC Pulm Med.* 2017 Dec 2;17(1):163.
 71. Zhang L, Xu Y, Li Y, Bao T, Gowd V, Chen W. Protective property of mulberry digest against oxidative stress-A potential approach to ameliorate dietary acrylamide-induced cytotoxicity. *Food Chem.* 2017 Sep 1;230:306-315.
 72. Feng M, Yin H, Peng H, Liu Z, Lu G, Dang Z. Hexavalent chromium induced oxidative stress and apoptosis in *Pycnoporus sanguineus*. *Environ Pollut.* 2017 Sep;228:128-139.
 73. Meng J, Lv Z, Qiao X, Li X, Li Y, Zhang Y, Chen C. The decay of Redox-stress Response Capacity is a substantive characteristic of aging: Revising the redox theory of aging. *Redox Biol.* 2017 Apr;11:365-374.
 74. Xuanfei L, Hao C, Zhujun Y, Yanming L, Jianping G. Imidazole 12 receptor inhibitor idazoxan regulates the progression of hepatic fibrosis via Akt-Nrf2-Smad2/3 signaling pathway. *Oncotarget.* 2017 Mar 28;8(13):21015-21030.
 75. Wang M, Wang L, Jia Z, Wang X, Yi Q, Zhao L, Song L. The versatile functions of LRR-only proteins in mollusk *Chlamys farreri*. *Dev Comp Immunol.* 2017 Dec;77:188-199.
 76. Hou B, Qiang G, Zhao Y, Yang X, Chen X, Yan Y, Wang X, Liu C, Zhang L, Du G. Salvianolic Acid A Protects Against Diabetic Nephropathy through Ameliorating Glomerular Endothelial Dysfunction via Inhibiting AGE-RAGE Signaling. *Cell Physiol Biochem.* 2017;44(6):2378-2394.
 77. Han X, Xue X, Zhao Y, Li Y, Liu W, Zhang J, Fan S. Rutin-Enriched Extract from *Coriandrum sativum* L. Ameliorates Ionizing Radiation-Induced Hematopoietic Injury. *Int J Mol Sci.* 2017 Apr 29;18(5). pii: E942.
 78. Wan B, Zhang Q, Ni J, Li S, Wen D, Li J, Xiao H, He P, Ou HY, Tao J, Teng Q, Lu J, Wu W, Yao YF. Type VI secretion system contributes to Enterohemorrhagic *Escherichia coli* virulence by secreting catalase against host reactive oxygen species (ROS). *PLoS Pathog.* 2017 Mar 13;13(3):e1006246.
 79. Lu QB, Wan MY, Wang PY, Zhang CX, Xu DY, Liao X, Sun HJ. Chicoric acid prevents PDGF-BB-induced VSMC dedifferentiation, proliferation and migration by suppressing ROS/NF κ B/mTOR/P70S6K signaling cascade. *Redox Biol.* 2018 Apr;14:656-668.
 80. Wang N, Zhang J, Qin M, Yi W, Yu S, Chen Y, Guan J, Zhang R. Amelioration of streptozotocin induced pancreatic β cell damage by morin: Involvement of the AMPK FOXO3 catalase signaling pathway. *Int J Mol Med.* 2018 Mar;41(3):1409-1418.
 81. Zhao L, Qi L, Li C, Li L, Jin L, Yuan J. SVCV impairs mitochondria complex III resulting in accumulation of hydrogen peroxide. *Fish Shellfish Immunol.* 2018 Apr;75:58-65.
 82. Feng M, Yin H, Cao Y, Peng H, Lu G, Liu Z, Dang Z. Cadmium-induced stress response of *Phanerochaete chrysosporium* during the biodegradation of 2,2',4,4'-tetrabromodiphenyl ether (BDE-47). *Ecotoxicol Environ Saf.* 2018 Jun 15;154:45-51.
 83. Tao J, Xu J, Chen F, Xu B, Gao J, Hu Y. Folate acid-Cyclodextrin/Docetaxel induces apoptosis in KB cells via the intrinsic mitochondrial pathway and displays antitumor activity in vivo. *Eur J Pharm Sci.* 2018 Jan 1;111:540-548.
 84. Wang J, Zhang J, Chen X, Yang Y, Wang F, Li W, Awuti M, Sun Y, Lian C, Li Z, Wang M, Xu JY, Jin C, Tian H, Gao F, Zhang J, Sinha D, Lu L, Xu GT. miR-365 promotes diabetic retinopathy through inhibiting Timp3 and increasing oxidative stress. *Exp Eye Res.* 2018 Mar;168:89-99.
 85. Zhang S, Huang Q, Wang Q, Wang Q, Cao X, Zhao L, Xu N, Zhuge Z, Mao J, Fu X, Liu R, Wilcox CS, Patzak A, Li L, Lai EY. Enhanced Renal Afferent Arteriolar Reactive Oxygen Species and Contractility to Endothelin-1 Are Associated with Canonical Wnt Signaling in Diabetic Mice. *Kidney Blood Press R.* 2018;43(3):860-871.
 86. Zhao L, Qi L, Li C, Li L, Jin L, Yuan J. SVCV impairs mitochondria complex III resulting in accumulation of hydrogen peroxide. *Fish Shellfish Immunol.* 2018 Apr;75:58-65.
 87. Lu QB, Wan MY, Wang PY, Zhang CX, Xu DY, Liao X, Sun HJ. Chicoric acid prevents PDGF-BB-induced VSMC dedifferentiation, proliferation and migration by suppressing ROS/NF κ B/mTOR/P70S6K signaling cascade. *Redox Biol.* 2018 Apr;14:656-668.
 88. Ma TJ, Lan DH, He SZ, Ye Z, Li P, Zhai W, Chen WQ, Huang Y, Fu Y, Sun A, Wang YB, Ye Z, Li JL, Gao Y, Yan XL, Li ZH. Nrf2 protects human lens epithelial cells against H₂O₂-induced oxidative and ER stress: The ATF4 may be involved. *Exp Eye Res.* 2018 Apr;169:28-37.
 89. Xue J, Gu H, Liu D, Ma W, Wei X, Zhao L, Liu Y, Zhang C, Yuan Z. Mitochondrial dysfunction is implicated in retinoic acid-induced spina bifida aperta in rat fetuses. *Int J Dev Neurosci.* 2018 Aug;68:39-44.
 90. Chen L, Zhang J, Li C, Wang Z, Li J, Zhao D, Wang S, Zhang H, Huang Y, Guo X. Glycine Transporter-1 and glycine receptor mediate the antioxidant effect of glycine in diabetic rat islets and INS-1 cells. *Free Radical Bio Med.* 2018 Aug 1;123:53-61.
 91. Wang XX, Zhang TY, Dao GH, Hu HY. Interaction between 1,2-benzisothiazol-3(2H)-one and microalgae: Growth inhibition and detoxification mechanism. *Aquat Toxicol.* 2018 Dec;205:66-75.
 92. Xu S, Zhao X, Liu F, Cao Y, Wang B, Wang X, Yin M, Wang Q, Feng

- XCrucial role of oxidative stress in bactericidal effect of parthenolide against *Xanthomonas oryzae* pv. *oryzae*. *Pest Manag Sci* . 2018 Dec;74(12):2716-2723.
93. Wu Z, Wang Y, Meng X, Wang X, Li Z, Qian S, Wei Y, Shu L, Ding Y, Wang P, Peng Y Total C-21 steroidal glycosides, isolated from the root tuber of *Cynanchum auriculatum* Royle ex Wight, attenuate hydrogen peroxide-induced oxidative injury and inflammation in L02 cells. *Int J Mol Med* . 2018 Dec;42(6):3157-3170
94. Liang Z, Yuan Z, Li G, Fu F, Shan Y Hypolipidemic, Antioxidant, and Antiapoptotic Effects of Polysaccharides Extracted from Reishi Mushroom, *Ganoderma lucidum* (Leysser: Fr) Karst, in Mice Fed a High-Fat Diet. *J Med Food* . 2018 Dec;21(12):1218-1227..
95. Wang Z, Zhou Y, Yu Y, He K, Cheng LM. Lipopolysaccharide preconditioning increased the level of regulatory B cells in the spleen after acute ischaemia/reperfusion in mice. *Brain Res* . 2018 Dec 15;1701:46-57.
96. Li Y, Li P, Wang T, Zhang FJ, Huang XX, Hou BK. The maize secondary metabolism glycosyltransferase UFGT2 modifies flavonols and contributes to plant acclimation to abiotic stresses. *Ann Bot-London* . 2018 Dec 31;122(7):1203-1217.
97. Ren C, Wu H, Li D, Yang Y, Gao Y, Jizhang Y, Liu D, Ji X, Zhang X Remote Ischemic Conditioning Protects Diabetic Retinopathy in Streptozotocin-induced Diabetic Rats via Anti-Inflammation and Antioxidation. *Aging Dis* . 2018 Dec 4;9(6):1122-1133.
98. Wang L, Wang F, Liu S, Yang X, Yang J, Ming D. VEGF attenuates 2-VO induced cognitive impairment and neuronal injury associated with the activation of PI3K/Akt and Notch1 pathway. *Exp Gerontol* . 2018 Feb;102:93-100.
99. Dai J, Gu L, Su Y, Wang Q, Zhao Y, Chen X, Deng H, Li W, Wang G, Li K. Inhibition of curcumin on influenza A virus infection and influenzal pneumonia via oxidative stress, TLR2/4, p38/JNK MAPK and NF- κ B pathways. *Int Immunopharmacol* . 2018 Jan;54:177-187.

Version 2020.03.20