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- umbilical vein endothelial cells against lysophosphatidylcholine-induced apoptosis by upregulating superoxide dismutase and glutathione peroxidase. *IUBMB Life* . 2014 Oct;66(10):711-22.
2. Zhao H, Wang R, Tao Z, Yan F, Gao L, Liu X, Wang N, Min L, Jia Y, Zhao Y, Ji X, Luo Y. Activation of T-LAK-cell-originated protein kinase-mediated antioxidation protects against focal cerebral ischemia-reperfusion injury. *FEBS J* . 2014 Oct;281(19):4411-20.
 3. Li X, Xu L, Zhou W, Zhao Q, Wang Y. Chronic exposure to microcystin-LR affected mitochondrial DNA maintenance and caused pathological changes of lung tissue in mice. *Environ Pollut* . 2015 Dec 16;210:48-56.
 4. Liu J, Tang X, Wang Y, Zang Y, Zhou B. A *Zostera marina* manganese superoxide dismutase gene involved in the responses to temperature stress. *Gene* . 2016 Jan 10;575(2 Pt 3):718-24.
 5. Chen W, Su H, Xu Y, Bao T, Zheng X. Protective effect of wild raspberry (*Rubus hirsutus* Thunb.) extract against acrylamide-induced oxidative damage is potentiated after simulated gastrointestinal digestion. *Food Chem* . 2016 Apr 1;196:943-52.
 6. Sun Y, Xiu C, Liu W, Tao Y, Wang J, Qu YL. Grape seed proanthocyanidin extract protects the retina against early diabetic injury by activating the Nrf2 pathway. *Exp Ther Med* . 2016 Apr;11(4):1253-1258.
 7. Lin J, Zou Y, Cao K, Ma C, Chen Z. The impact of heterologous catalase expression and superoxide dismutase overexpression on enhancing the oxidative resistance in *Lactobacillus casei*. *J IND MICROBIOL BIOT* . 2016 May;43(5):703-11.
 8. Wang J, Chen X, Wang F, Zhang J, Li P, Li Z, Xu J, Gao F, Jin C, Tian H, Zhang J, Li W, Lu L, Xu GT. OFD1, as a Ciliary Protein, Exhibits Neuroprotective Function in Photoreceptor Degeneration Models. *PLoS One* . 2016 May 19;11(5):e0155860.
 9. Zhuang C, Xu NW, Gao GM, Ni S, Miao KS, Li CK, Wang LM, Xie HG. Polysaccharide from *Angelica sinensis* protects chondrocytes from H₂O₂-induced apoptosis through its antioxidant effects in vitro. *Int J Biol Macromol* . 2016 Jun;87:322-8.
 10. Guo H, Zhang N, Liu D, Wang P, Ma X. Inhibitory effect on the proliferation of human hepatoma induced by cell-permeable Biomed Pharmacother . 2016 Oct;83:1379-1386.
 11. Yang J, Liu X, Zhang X, Jin Q, Li J. Phenolic Profiles, Antioxidant Activities, and Neuroprotective Properties of Mulberry. *J Food Sci* . 2016 Oct;81(10):C2439-C2446.
 12. Sun X, Jiao X, Ma Y, Liu Y, Zhang L, He Y, Chen Y. Trimethylamine N-oxide induces inflammation and endothelial dysfunction in human umbilical vein endothelial cells via activating ROS-TXNIP-NLRP3 inflammasome. *BIOCHEM BIOPH RES CO* . 2016 Dec 2;481(1-2):63-70.
 13. Pan K, Li X, Chen Y, Zhu D, Li Y, Tao G, Zuo Z. Deferoxamine pre-treatment protects against postoperative cognitive dysfunction of aged rats by depressing microglial activation via ameliorating iron accumulation in hippocampus. *Neuropharmacology* . 2016 Dec;111:180-194.
 14. Dong K, Hao P, Xu S, Liu S, Zhou W, Yue X, Rausch-Fan X, Liu Z. Alpha-Lipoic Acid Alleviates High-Glucose Suppressed Osteogenic Differentiation of MC3T3-E1 Cells via Antioxidant Effect and PI3K/Akt Signaling Pathway. *CELL PHYSIOL BIOCHEM* . 2017;42(5):1897-1906.
 15. Zhang W, Zhang Y, Guo X, Zeng Z, Wu J, Liu Y, He J, Wang R, Huang Q, Chen Z. Sirt1 Protects Endothelial Cells against LPS-Induced Barrier Dysfunction. *Oxid Med Cell Longev* . 2017;2017:4082102.
 16. Cao XN, Shen LJ, Wu SD, Yan C, Zhou Y, Xiong G, Wang YC, Liu Y, Liu B, Tang XL, Guo M, Liu DY, Long CL, Sun M, He DW, Lin T, Wei GH. Urban fine particulate matter exposure causes male reproductive injury through destroying blood-testis barrier (BTB) integrity. *Toxicol Lett* . 2017 Jan 15;266:1-12.
 17. Huang W, Guo HL, Deng X, Zhu TT, Xiong JF, Xu YH, Xu Y. Short-Chain Fatty Acids Inhibit Oxidative Stress and Inflammation in Mesangial Cells Induced by High Glucose and Lipopolysaccharide. *EXP CLIN ENDOCR DIAB* . 2017 Jan 3. doi: 10.1055/s-0042-121493.
 18. Chen W, Su H, Xu Y, Jin C. In vitro gastrointestinal digestion promotes the protective effect of blackberry extract against acrylamide-induced oxidative stress. *SCI REP-UK* . 2017 Jan 13;7:40514.
 19. Cao XN, Shen LJ, Wu SD, Yan C, Zhou Y, Xiong G, Wang YC, Liu Y, Liu B, Tang XL, Guo M, Liu DY, Long CL, Sun M, He DW, Lin T, Wei GH. Urban fine particulate matter exposure causes male reproductive injury through destroying blood-testis barrier (BTB) integrity. *Toxicol Lett* . 2017 Jan 15;266:1-12.
 20. Hu W, Wang H, Liu Z, Liu Y, Wang R, Luo X, Huang Y. Neuroprotective effects of lycopene in spinal cord injury in rats via antioxidative and anti-apoptotic pathway. *Neurosci Lett* . 2017 Feb 3;642:107-112.
 21. Li H, Li Z, Peng L, Jiang N, Liu Q, Zhang E, Liang B, Li R, Zhu H. Lycium barbarum polysaccharide protects human keratinocytes against UVB-induced photo-damage. *FREE RADICAL RES* . 2017 Feb;51(2):200-210.
 22. Hu W, Wang H, Liu Z, Liu Y, Wang R, Luo X, Huang Y. Neuroprotective effects of lycopene in spinal cord injury in rats via antioxidative and anti-apoptotic pathway. *Neurosci Lett* . 2017 Mar 6;642:107-112.
 23. Yuan J, Zhu C, Hong Y, Sun Z, Fang X, Wu B, Li S. The role of cPLA2 in Methylglyoxal-induced cell apoptosis of HUVECs. *TOXICOL APPL PHARM* . 2017 May 15;323:44-52.
 24. Gong W, Li J, Chen Z, Huang J, Chen Q, Cai W, Liu P, Huang H. Polydatin promotes Nrf2-ARE anti-oxidative pathway through activating CKIP-1 to resist HG-induced up-regulation of FN and ICAM-1 in GMCs and diabetic mice kidneys. *FREE RADICAL BIO MED* . 2017 May;106:393-405.
 25. Wang M, Zhang Y, Guo P. Effect of florfenicol and thiamphenicol exposure on the photosynthesis and antioxidant system of *Microcystis flos-aquae*. *Aquat Toxicol* . 2017 May;186:67-76.
 26. Chang Y, Li S, Guo W, Yang Y, Zhang W, Zhang Q, He Y, Yi X, Cui T, An Y, Song P, Jian Z, Liu L, Li K, Wang G, Gao T, Wang L, Li C. Simvastatin Protects Human Melanocytes from H₂O₂-Induced Oxidative Stress by Activating Nrf2. *J Invest Dermatol* . 2017 Jun;137(6):1286-1296.
 27. Jiao YH, Dou M, Wang G, Li HY, Liu JS, Yang X, Yang WD. Exposure of okadaic acid alters the angiogenesis in developing chick embryos. *Toxicol* . 2017 Jul;133:74-81.
 28. Leng ZG, Lin SJ, Wu ZR, Guo YH, Cai L, Shang HB, Tang H, Xue YJ, Lou MQ, Zhao W, Le WD, Zhao WG, Zhang X, Wu ZB. Activation of DRD5 (dopamine receptor D5) inhibits tumor growth by autophagic cell death. *Autophagy* . 2017 Aug 3;13(8):1404-1419.
 29. Fu Q, Huang T, Wang X, Lu C, Liu F, Yang G, Wang Y, Wang B. Association of elevated reactive oxygen species and hyperthermia induced radiosensitivity in cancer stem-like cells. *ONCOTARGET* . 2017 Oct 9;8(60):101560-101571.
 30. Peng J, Deng X, Huang W, Yu JH, Wang JX, Wang JP, Yang SB, Liu X, Wang L, Zhang Y, Zhou XY, Yang H, He YZ, Xu FY. Irisin protects against neuronal injury induced by oxygen-glucose deprivation in part depends on the inhibition of ROS-NLRP3 inflammatory signaling pathway. *Mol Immunol* . 2017 Nov;91:185-194.
 31. Dong X, Niu Y, Ding Y, Wang Y, Zhao J, Leng W, Qin L. Formulation and Drug Loading Features of Nano-Erythrocytes. *Nanoscale Res Lett* . 2017 Dec;12(1):202.
 32. Liu SY, Lu S, Yu XL, Yang SG, Liu W, Liu XM, Wang SW, Zhu J, Ji M, Liu DQ, Zhang ZP, Liu RT. Fruitless Wolfberry-Sprout Extract Rescued Cognitive Deficits and Attenuated Neuropathology in Alzheimer's Disease Transgenic Mice. *Curr Alzheimer Res* . 2018;15(9):856-868.
 33. Lu MC, Jiao Q, Liu T, Tan SJ, Zhou HS, You QD, Jiang ZY. Discovery of a head-to-tail cyclic peptide as the Keap1-Nrf2 protein-protein interaction inhibitor with high cell potency. *Eur J Med Chem* . 2018 Jan 1;143:1578-1589.
 34. Fu Q, Li C, Yu L. Gambogic acid inhibits spinal cord injury and inflammation through suppressing the p38 and Akt signaling pathways. *Mol Med Rep* . 2018 Jan;17(1):2026-2032.
 35. Wang M, Wang B, Jiang K, Liu M, Shi X, Wang L. A mitochondrial manganese superoxide dismutase involved in innate immunity is essential for the survival of *Chlamys farreri*. *FISH SHELLFISH IMMUN* . 2018 Jan;72:282-290.
 36. Ge J, Chen L, Yang Y, Lu X, Xiang Z. Sparstolonin B prevents lumbar intervertebral disc degeneration through toll like receptor 4, NADPH oxidase activation and the protein kinase B signaling pathway. *Mol Med Rep* . 2018 Jan;17(1):1347-1353.
 37. Jiang Y, Wang W, Liu ZY, Xie Y, Qian Y, Cai XN. Overexpression of miR-130a-3p/301a-3p attenuates high glucose-induced MPC5 podocyte dysfunction through suppression of TNF- α signaling. *Exp Ther Med* . 2018 Jan;15(1):1021-1028.
 38. Dai J, Gu L, Su Y, Wang Q, Zhao Y, Chen X, Deng H, Li W, Wang G, Li K. Inhibition of curcumin on influenza A virus infection and influenzal pneumonia via oxidative stress, TLR2/4, p38/JNK MAPK and NF- κ B pathways. *Int Immunopharmacol* . 2018 Jan;54:177-187.

39. Nie HT,Guo YX,Yao XL,Ma TW,Deng KP,Wang Z,Zhang GM,Sun LW,Wang ZY,Wang HC,Wang F. Acute nutrient treatment causes alterations in intra-follicular antioxidation and AKT signaling. *Reproduction* . 2018 Jan;155(1):37-49.
40. Gong W, Chen Z, Zou Y, Zhang L, Huang J, Liu P, Huang H. CKIP-1 affects the polyubiquitination of Nrf2 and Keap1 via mediating Smurf1 to resist HG-induced renal fibrosis in GMCs and diabetic mice kidneys. *FREE RADICAL BIO MED* . 2018 Feb 1;115:338-350.
41. Wang L,Wang F,Liu S,Yang X,Yang J,Ming D. VEGF attenuates 2-VO induced cognitive impairment and neuronal injury associated with the activation of PI3K/Akt and Notch1 pathway. *Exp Gerontol* . 2018 Feb;102:93-100.
42. Chen L,Yao H,Chen X,Wang Z,Xiang Y,Xia J,Liu Y,Wang Y. Ginsenoside Rg1 Decreases Oxidative Stress and Down-Regulates Akt/mTOR Signalling to Attenuate Cognitive Impairment in Mice and Senescence of Neural Stem Cells Induced by D-Galactose. *Neurochem Res* . 2018 Feb;43(2):430-440.
43. Ma J,Song D,Zhang Y,Chen L,Zhang S,Jia J,Chen T,Guo C,Tian L,Gao A,Niu P. SIRT1 exhibits antioxidative effects in HT22 cells induced by tert-butyl alcohol. *Environ Toxicol* . 2018 Feb;33(2):142-148.
44. Wang J, Zhang J, Chen X, Yang Y, Wang F, Li W, Awuti M, Sun Y, Lian C, Li Z, Wang M, Xu JY, Jin C, Tian H, Gao F, Zhang J, Sinha D, Lu L, Xu GT. miR-365 promotes diabetic retinopathy through inhibiting Timp3 and increasing oxidative stress. *Exp Eye Res* . 2018 Mar;168:89-99.
45. Zhang N,Feng H,Liao HH,Chen S,Yang Z,Deng W,Tang QZ. Myricetin attenuated LPS induced cardiac injury in vivo and in vitro. *Phytother Res* . 2018 Mar;32(3):459-470.
46. Cui K,Tang Z,Li CC,Wang T,Rao K,Wang SG,Liu JH,Chen Z. Lipoxin A4 improves erectile dysfunction in rats with type I diabetes by inhibiting oxidative stress and corporal fibrosis. *Asian J Androl* . 2018 Mar-Apr;20(2):166-172.
47. Xie X,Chen Q,Tao J. Astaxanthin Promotes Nrf2/ARE Signaling to Inhibit HG-Induced Renal Fibrosis in GMCs. *Mar Drugs* . 2018 Apr 5;16(4). pii: E117.
48. Yang J,Zhang Y,Zang G,Wang T,Yu Z,Wang S,Tang Z,Liu J. Adipose-derived stem cells improve erectile function partially through the secretion of IGF-1, bFGF, and VEGF in aged rats. *Andrology* . 2018 May;6(3):498-509.
49. Fang Z,Luo W,Luo Y. Protective effect of α -mangostin against CoCl₂-induced apoptosis by suppressing oxidative stress in H9C2 rat cardiomyoblasts. *Mol Med Rep* . 2018 May;17(5):6697-6704.
50. Wu D,Han R,Deng S,Liu T,Zhang T,Xie H,Xu Y. Protective Effects of Flagellin A N/C Against Radiation-Induced NLR Pyrin Domain Containing 3 Inflammasome-Dependent Pyroptosis in Intestinal Cells. *INT J RADIAT ONCOL* . 2018 May 1;101(1):107-117.
51. Wang X,Chen B,Sun J,Jiang Y,Zhang H,Zhang P,Fei B,Xu Y. Iron-induced oxidative stress stimulates osteoclast differentiation via NF- κ B signaling pathway in mouse model. *Metabolism* . 2018 Jun;83:167-176.
52. Chen M,Zheng J,Liu G,Xu E,Wang J,Fuqua BK,Vulpe CD,Anderson GJ,Chen H. Ceruloplasmin and hephaestin jointly protect the exocrine pancreas against oxidative damage by facilitating iron efflux. *Redox Biol* . 2018 Jul;17:432-439.
53. Pan L,Zhou L,Yin W,Bai J,Liu R. miR-125a induces apoptosis, metabolism disorder and migration impairment in pancreatic cancer cells by targeting Mfn2-related mitochondrial fission. *Int J Oncol* . 2018 Jul;53(1):124-136.
54. Zhang X,Li J,Cheng Y,Yi J,Liu X,Cheng W. Downregulation of CUEDC2 prevents doxorubicin-induced cardiotoxicity in H9c2 cells. *Mol Med Rep* . 2018 Jul;18(1):855-863.
55. Wang Y,Chen G,Yan J,Chen X,He F,Zhu C,Zhang J,Lin J,Pan G,Yu J,Pei M,Yang H,Liu T. Upregulation of SIRT1 by Kartogenin Enhances Antioxidant Functions and Promotes Osteogenesis in Human Mesenchymal Stem Cells. *Oxid Med Cell Longev* . 2018 Jul 15;2018:1368142.
56. Ding W,Zhao Y,Xu JW,Zhao P,Li T,Ma H,Reiter RJ,Yu X. Melatonin: A Multifunctional Molecule That Triggers Defense Responses against High Light and Nitrogen Starvation Stress in *Haematococcus pluvialis*. *J AGR FOOD CHEM* . 2018 Jul 25;66(29):7701-7711.
57. Yanfei Cui,Shengya Yang. Overexpression of Annexin A1 Protects Against Benzof[a]pyrene-induced Bronchial Epithelium Injury *Mol Med Rep*. 2018 Jul;18(1):349-357. doi: 10.3892/mmr.2018.8998.
58. Jin Y,Tu Q,Liu M. MicroRNA-125b regulates Alzheimer's disease through SphK1 regulation. *Mol Med Rep* . 2018 Aug;18(2):2373-2380.
59. Ma X,Hu B,Zou C,Han A,Xu Z,Zhang T,Yu W. The effects of hyperoxia liquid regulate cardiopulmonary bypass-induced myocardial damage through the Nrf2-ARE signaling pathway. *Mol Med Rep* . 2018 Aug;18(2):2342-2348.
60. Chen L,Zhang J,Li C,Wang Z,Li J,Zhao D,Wang S,Zhang H,Huang Y,Guo X. Glycine Transporter-1 and glycine receptor mediate the antioxidant effect of glycine in diabetic rat islets and INS-1 cells. *FREE RADICAL BIO MED* . 2018 Aug 1;123:53-61.
61. Hu Y,Pan J,Xin Y,Mi X,Wang J,Gao Q,Luo H. Gene Expression Analysis Reveals Novel Gene Signatures Between Young and Old Adults in Human Prefrontal Cortex. *Front Aging Neurosci* . 2018 Aug 27;10:259.
62. Liu H,Zhang X,Zhang S,Huang H,Wu J,Wang Y,Yuan L,Liu C,Zeng X,Cheng X,Zhuang D,Zhang H. Oxidative Stress Mediates Microcystin-LR-Induced Endoplasmic Reticulum Stress and Autophagy in KK-1 Cells and C57BL/6 Mice Ovaries. *Front Physiol* . 2018 Aug 6;9:1058.
63. Huang Y,Hu Z. UBIAD1 protects against oxygen-glucose deprivation/reperfusion-induced multiple subcellular organelles injury through PI3K/AKT pathway in N2A cells. *J Cell Physiol* . 2018 Sep;233(9):7480-7496.
64. Bao J,Ye C,Zheng Z,Zhou Z. Fmr1 protects cardiomyocytes against lipopolysaccharide-induced myocardial injury. *Exp Ther Med* . 2018 Sep;16(3):1825-1833.
65. Wang L,Chen HC,Yang X,Tao JJ,Liang G,Wu JZ,Wu WC,Wang Y,Song ZM,Zhang X. The novel chalcone analog L2H17 protects retinal ganglion cells from oxidative stress-induced apoptosis. *Neural Regen Res* . 2018 Sep;13(9):1665-1672.
66. Xu E,Chen M,Zheng J,Maimaitiming Z,Zhong T,Chen H. Deletion of hephaestin and ceruloplasmin induces a serious systemic iron deficiency and disrupts iron homeostasis. *BIOCHEM BIOPH RES CO* . 2018 Sep 10;503(3):1905-1910.
67. Ji C,Luo Y,Zou C,Huang L,Tian R,Lu Z. Effect of astragaloside IV on indoxyl sulfate-induced kidney injury in mice via attenuation of oxidative stress. *BMC PHARMACOL TOXICO* . 2018 Sep 3;19(1):53.
68. Fu Y,Wu HQ,Cui HL,Li YY,Li CZ. Gastroprotective and anti-ulcer effects of oxymatrine against several gastric ulcer models in rats: Possible roles of antioxidant, antiinflammatory, and prosurvival mechanisms. *Phytother Res* . 2018 Oct;32(10):2047-2058.
69. Liu S,Yuan J,Yue W,Bi Y,Shen X,Gao J,Xu X,Lu Z. GCN2 deficiency protects against high fat diet induced hepatic steatosis and insulin resistance in mice. *BBA-MOL BASIS DIS* . 2018 Oct;1864(10):3257-3267.
70. Jiang H,Gao X,Gong J,Yang Q,Lan R,Wang T,Liu J,Yin C,Wang S,Liu Z. Downregulated Expression of Solute Carrier Family 26 Member 6 in NRK-52E Cells Attenuates Oxalate-Induced Intracellular Oxidative Stress. *Oxid Med Cell Longev* . 2018 Oct 10;2018:1724648.
71. Lu C,Chen X,Wang Q,Xu X,Xu B. TNF α promotes glioblastoma A172 cell mitochondrial apoptosis via augmenting mitochondrial fission and repression of MAPK-ERK-YAP signaling pathways. *ONCOTARGETS THER* . 2018 Oct 18;11:7213-7227.
72. Liu Q,Hu Y,Zhang M,Yan Y,Yu H,Ge L. microRNA-451 protects neurons against ischemia/reperfusion injury-induced cell death by targeting CELF2. *Neuropsychiatr Dis Treat* . 2018 Oct 23;14:2773-2782.
73. Deng J,Liu Q,Zhang Q,Zhang C,Liu D,Fan D,Yang H. Comparative study on composition, physicochemical and antioxidant characteristics of different varieties of kiwifruit seed oil in China. *Food Chem* . 2018 Oct 30;264:411-418.
74. Zheng JW,Liu SL,Lu SH,Li HY,Liu JS,Yang WD. Proteomic profile in the mussel *Perna viridis* after short-term exposure to the brown tide alga *Aureococcus anophagefferens*. *ECOTOX ENVIRON SAFE* . 2018 Oct 30;162:365-375.
75. Xu N,Kan P,Yao X,Yang P,Wang J,Xiang L,Zhu Y. Astragaloside IV reversed the autophagy and oxidative stress induced by the intestinal microbiota of AIS in mice. *J Microbiol* . 2018 Nov;56(11):838-846.
76. Zhang D,Li Y,Zhang T,Liu J,Jahejo AR,Yang L,Chen P,Ning G,Huo N,Ma H,Yan F,Tian W. Protective effects of zinc and N-acetyl-L-cysteine supplementation against cadmium induced erythrocyte cytotoxicity in Arbor Acres broiler chickens (*Gallus gallus domesticus*). *ECOTOX ENVIRON SAFE* . 2018 Nov 15;163:331-339.
77. Yao S,Yan W. Overexpression of Mst1 reduces gastric cancer cell viability by repressing the AMPK-Sirt3 pathway and activating mitochondrial fission. *ONCOTARGETS THER* . 2018 Nov

- 29;11:8465-8479.
78. Chen Q, Tao J, Li G, Zheng D, Tan Y, Li R, Tian L, Li Z, Cheng H, Xie X. Astaxanthin ameliorates experimental diabetes-induced renal oxidative stress and fibronectin by upregulating connexin43 in glomerular mesangial cells and diabetic mice. *Eur J Pharmacol*. 2018 Dec 5;840:33-43.
 79. Lin S, Ren A, Wang L, Santos C, Huang Y, Jin L, Li Z, Greene NDE. Aberrant methylation of Pax3 gene and neural tube defects in association with exposure to polycyclic aromatic hydrocarbons. *Clin Epigenetics*. 2019 Jan 21;11(1):13.
 80. Yang J, Zhou X, Zeng X, Hu O, Yi L, Mi M. Resveratrol attenuates oxidative injury in human umbilical vein endothelial cells through regulating mitochondrial fusion via TyrRS-PARP1 pathway. *Nutr Metab (Lond)*. 2019 Jan 30;16:9.
 81. Zhang H, Chen Y, Pei Z, Gao H, Shi W, Sun M, Xu Q, Zhao J, Meng W, Xiao K. Protective effects of polydatin against sulfur mustard-induced hepatic injury. *Toxicol Appl Pharmacol*. 2019 Mar 15;367:1-11.
 82. Wang SQ, Yang XY, Cui SX, Gao ZH, Qu XJ. Heterozygous knockout insulin-like growth factor-1 receptor (IGF-1R) regulates mitochondrial functions and prevents colitis and colorectal cancer. *Free Radic Biol Med*. 2019 Apr;134:87-98.
 83. Chang G, Chen Y, Zhang H, Zhou W. Trans sodium crocetininate alleviates ischemia/reperfusion-induced myocardial oxidative stress and apoptosis via the SIRT3/FOXO3a/SOD2 signaling pathway. *Int Immunopharmacol*. 2019 Jun;71:361-371.
 84. Xu X, Huang L, Zhang Z, Tong J, Mi J, Wu Y, Zhang C, Yan H. Targeting non-oncogene ROS pathway by alantolactone in B cell acute lymphoblastic leukemia cells. *Life Sci*. 2019 Jun 15;227:153-165.
 85. Rao G, Zhang W, Song S. MicroRNA-217 inhibition relieves cerebral ischemia/reperfusion injury by targeting SIRT1. *Mol Med Rep*. 2019 Aug;20(2):1221-1229.
 86. Li X, Qi M, Sun X, Weir MD, Tay FR, Oates TW, Dong B, Zhou Y, Wang L, Xu HHK. Surface treatments on titanium implants via nanostructured ceria for antibacterial and anti-inflammatory capabilities. *Acta Biomater*. 2019 Aug;94:627-643.
 87. Li Y, Ma S, Zhang Y, Yao M, Zhu X, Guan F. (-)-Epicatechin mitigates radiation-induced intestinal injury and promotes intestinal regeneration via suppressing oxidative stress. *Free Radic Res*. 2019 Aug;53(8):851-864.
 88. Lu MC, Zhao J, Liu YT, Liu T, Tao MM, You QD, Jiang ZY. CPUY192018, a potent inhibitor of the Keap1-Nrf2 protein-protein interaction, alleviates renal inflammation in mice by restricting oxidative stress and NF- κ B activation. *Redox Biol*. 2019 Sep;26:101266.
 89. Liu C, Wu QQ, Cai ZL, Xie SY, Duan MX, Xie QW, Yuan Y, Deng W, Tang QZ. Zingerone attenuates aortic banding-induced cardiac remodeling via activating the eNOS/Nrf2 pathway. *J Cell Mol Med*. 2019 Sep;23(9):6466-6478.
 90. Zhang J, Zhang SD, Wang P, Guo N, Wang W, Yao LP, Yang Q, Efferth T, Jiao J, Fu YJ. Pinolenic acid ameliorates oleic acid-induced lipogenesis and oxidative stress via AMPK/SIRT1 signaling pathway in HepG2 cells. *Eur J Pharmacol*. 2019 Oct 15;861:172618.
 91. Song Y, Li X, Gong X, Zhao X, Ma Z, Xia T, Gu X. Green tea polyphenols improve isoflurane-induced cognitive impairment via modulating oxidative stress. *J Nutr Biochem*. 2019 Nov;73:108213.
 92. Chen J, Huang Z, Wu X, Kang J, Ren Y, Gao W, Lu X, Wang J, Ding W, Nakabeppu Y, Fan Y, Wang Y. Oxidative stress induces different tissue dependent effects on Mutyh-deficient mice. *Free Radic Biol Med*. 2019 Nov 1;143:482-493.
 93. Li X, Qi M, Li C, Dong B, Wang J, Weir MD, Imazato S, Du L, Lynch CD, Xu L, Zhou Y, Wang L, Xu HHK. Novel nanoparticles of cerium-doped zeolitic imidazolate frameworks with dual benefits of antibacterial and anti-inflammatory functions against periodontitis. *J Mater Chem B*. 2019 Nov 28;7(44):6955-6971.
 94. Tang J, Xiong L, Shu X, Chen W, Li W, Li J, Ma L, Xiao Y, Li L. Antioxidant effects of bioactive compounds isolated from cordyceps and their protective effects against UVB-irradiated HaCaT cells. *J Cosmet Dermatol*. 2019 Dec;18(6):1899-1906.
 95. Jiang H, Li R, Zhang Z, Chang C, Liu Y, Liu Z, He Q, Wang Q. Retinoid X receptor α (RXR α)-mediated erythroid-2-related factor-2 (NRF2) inactivation contributes to N,N-dimethylformamide (DMF)-induced oxidative stress in HL-7702 and HuH6 cells. *J Appl Toxicol*. 2020 Apr;40(4):470-482.

Version 2021.01.21