



碧云天网站



微信公众号

碧云天生物技术/Beyotime Biotechnology

订货热线: 400-168-3301或800-8283301

订货e-mail: order@beyotime.com

技术咨询: info@beyotime.com

网址: http://www.beyotime.com

MTT

产品编号	产品名称	包装
ST316	MTT	500mg

产品简介:

- MTT即Methylthiazolyldiphenyl-tetrazolium bromide, 也称Thiazolyl blue tetrazolium bromide, 中文名为噻唑蓝。分子式为 $C_{18}H_{16}BrN_5S$, 分子量为414.32, CAS Number 298-93-1, Ultra Pure Grade。进口分装。
- 本产品为黄色至橙黄粉末。
- 常用于基于MTT检测方法的细胞增殖或细胞毒性检测。
- 溶于水或DMSO。常配制成5mg/ml的储备液。

包装清单:

产品编号	产品名称	包装
ST316	MTT	500mg
—	说明书	1份

保存条件:

4°C避光保存。

注意事项:

- 本产品仅限于专业人员的科学研究用, 不得用于临床诊断或治疗, 不得用于食品或药品, 不得存放于普通住宅内。
- 为了您的安全和健康, 请穿实验服并戴一次性手套操作。

使用本产品的文献:

1. Fan QY, Liu QH, Li CP, Wang J. Experimental study on the effect of proto porphyrin disodium in vitro on anti-hbv. Modern Preventive Medicine. 2007;34(19):3614-5.
2. Yang H, Chen J, Yang G, Zhang XH, Li Y, Wang M. Characterization and pathogenicity of the zinc metalloprotease empA of *Vibrio anguillarum* expressed in *Escherichia coli*. Curr Microbiol. 2007 Mar;54(3):244-8.
3. Li ZJ, Zhang K, Lou SQ, Zheng YF. A Cellular Study on Cytotoxicity of Magnesium-calcium Alloy. Chinese journal of bone and joint injury. 2007 Sep;22(9).
4. Xu J, Sun Y, Huang J, Chen C, Liu G, Jiang Y, Zhao Y, Jiang Z. Photo killing cancer cells using highly cell-specific antibody-TiO₂ bio conjugates and electroporation. Bio electro chemistry. 2007 Nov;71(2):217-22.
5. Zhan XD, Wang KX, Li CP. Studies on Anti-HBV Effect in vitro of Snail Polysaccharide. Chinese Journal of Experimental Traditional Medical Formulae. 2006 ;17(11):21912-2192.
6. Bai L, Pang WJ, Yang YJ, Yang GS. Modulation of Sirt1 by resveratrol and nicotinamide alters proliferation and differentiation of pig preadipocytes. Mol Cell Bio chem. 2008 Jan;307(1-2):129-40.
7. Zhou XM, Xu GX, Zhao DM. In vitro effect of prion peptide PrP¹⁰⁶⁻¹²⁶ on mouse macrophages: Possible role of macrophages in transport and proliferation for prion protein. Microb Pathog. 2008 Feb;44(2):129-34.
8. Tian Y, Wang Q, Liu Q, Ma Y, Cao X, Zhang Y. Role of RpoS in stress survival, synthesis of extracellular autoinducer 2, and virulence in *Vibrio alginolyticus*. Arch Microbiol. 2008 Nov;190(5):585-94.
9. Zhang SH, Wang WQ, Wang JL. Protective effect of tetrahydroxystilbene glucoside on cardiotoxicity induced by doxorubicin in vitro and in vivo. Acta Pharmacol Sin. 2009;30(11):1479-87.
10. Lu S, Xia D, Huang G, Jing H, Wang Y, Gu H. Concentration effect of gold nanoparticles on proliferation of keratinocytes. Colloids Surf B Biointerfaces. 2010;81(2):406-11.
11. Cui J, Ou S, He WJ, Du L, Zhao YD, Ruan HZ. Prevention of extracellular ADP-induced ATP accumulation of the cultured rat spinal astrocytes via P2Y₁-mediated inhibition of AMPK. Neurosci Lett. 2011 Oct 10;503(3):244-9.
12. Li S, Tang Y, Zhang J, Guo X, Shen H. 3A4, a new potential target for B and myeloid lineage leukemias. J Drug Target. 2011 Nov;19(9):797-804.
13. Xiao H, Lv F, Xu W, Zhang L, Jing P, Cao X. Deprenyl prevents MPP(+)-induced oxidative damage in PC12 cells by the upregulation of Nrf2-mediated NQO1 expression through the activation of PI3K/Akt and Erk. Toxicology. 2011 Dec 18;290(2-3):286-94.
14. Liao H, Qi R, Shen M, Cao X, Guo R, Zhang Y, Shi X. Improved cellular response on multiwalled carbon nanotube-incorporated electrospun polyvinylalcohol/chitosan nanofibrous scaffolds. Colloids Surf B Biointerfaces. 2011 Jun 1;84(2):528-35.
15. L Yu-liang, W Kai-bin, L Hong, H Yu-qin, Z Feng, L Xin-gang. Olfactory ensheathing cell viability in different cryopreservation systems* ★ Journal of Clinical Rehabilitative Tissue Engineering Research. 2011 Nov;15(46).
16. Lu J, Xu SY, Zhang QG, Lei HY. Bupivacaine induces reactive oxygen species production via activation of the AMP-activate dprotein kinase-dependent pathway. Pharmacology. 2011;87(3-4):121-9.
17. Chen QY, Liang ZQ, Sun T, Sun XJ. Grape skin polyphenols (GSPs) inhibit the migration of mouse mammary carcinoma 4T1 cells in Vitro. CLINICAL ONCOLOGY AND CANCER RESEARCH. 2011; 8(1):33-37.
18. Wang J, Sun P, Bao Y, Dou B, Song D, Li Y. Vitamin E renders protection to PC12 cells against oxidative damage and apoptosis induced by single-walled carbon nanotubes. Toxicol In Vitro. 2012 Feb; 26(1):32-41.
19. Cui Y, Zhao Y, Tian Y, Zhang W, Lü X, Jiang X. The molecular mechanism of action of bactericidal gold nanoparticles on *Escherichia coli*. Biomaterials. 2012 Mar;33(7):2327-33.
20. Sun LC, Lin JT, Li W, Zhang L, Zhou TL, Zhang XY. Nicotine inhibits histone deacetylase 6 activity and chaperone-dependent activation of the glucocorticoid receptor in A549 cells. Chin Med J (Engl). 2012 Feb; 125(4):662-6.

21. Liu M, Yang S, Jin L, Hu D, Wu Z, Yang S. Chemical constituents of the ethyl acetate extract of *Belamcanda chinensis* (L.) DC roots and their antitumor activities. *Molecules*. 2012 May 24;17(5):6156-69.
22. Sun T, Chen QY, Wu LJ, Yao XM, Sun XJ. Antitumor and anti metastatic activities of grape skin polyphenols in a murine model of breast cancer. *Food Chem Toxicol*. 2012 Oct;50(10):3462-7.
23. Lü L, Tang D, Wang L, Huang LQ, Jiang GS, Xiao XY, Zeng FQ. Gambogic acid inhibits TNF- α -induced invasion of human prostate cancer PC3 cells in vitro through PI3K/Akt and NF- κ B signaling pathways. *Acta Pharmcol Sin*. 2012 Apr;33(4):531-41.
24. Xu S, Dong D, Wang WJ, Fu L, Xu A. Transferable crosslinked chitosan membranes for human melanocyte culture. *J Biomed Mater Res A*. 2012 Mar;100(3):673-83.
25. Fang X, Wu P, Li J, Qi L, Tang Y, Jiang W, Zhao S. Combination of apoptin with photodynamic therapy induces nasopharyngeal carcinoma cell death in vitro and in vivo. *Oncol Rep*. 2012 Dec;28(6):2077-82.
26. Li Z, Li B, Wang M, Xie M, Shen H, Shen S, Wang X, Guo X, Yao M, Jin Y. The role of endosome evasion bypass in the reversal of multidrug resistance by lipid/nanoparticle assemblies. *J Mater Chem B*. 2012 Jul 31;10:1466-74.
27. Wen X, Xu S, Liu H, Zhang Q, Liang H, Yang C, Wang H. Neurotoxicity induced by bupivacaine via T-type calcium channels in SH-SY5Y cells. *PLoS One*. 2013 May 2;8(5):e62942.
28. Yang S, Zhao Q, Xiang H, Liu M, Zhang Q, Xue W, Song B, Yang S. Antiproliferative activity and apoptosis-inducing mechanism of constituents from *Toona sinensis* on human cancer cells. *Cancer Cell Int*. 2013 Feb 9;13(1):12.
29. Gao C, Wang X, Chen L, Wang JH, Gao ZT, Wang H. Knockdown of Bcl-3 inhibits cell growth and induces DNA damage in HTLV-1-infected cells. *Asian Pac J Cancer Prev*. 2013;14(1):405-8.
30. Gao Y, Liu Z, Zhang X, He J, Pan Y, Hao F, Xie L, Li Q, Qiu X, Wang E. Inhibition of cytoplasmic GSK-3 β increases cisplatin resistance through activation of Wnt/ β -catenin signaling in A549/DDP cells. *Cancer Lett*. 2013 Aug 9;336(1):231-9.
31. Xu Y, Wang N, Ding Y, Wang C, Yu Y, Liu S, Wang X, Li Z. Ammonium chloride enhances cisplatin cytotoxicity through DNA double-strand breaks in human cervical cancer cells. *Oncol Rep*. 2013 Sep;30(3):1195-200.
32. Li JZ, Wu JH, Yu SY, Shao QR, Dong XM. Inhibitory effects of paeoniflorin on lysophosphatidylcholine-induced inflammatory factor production in human umbilical vein endothelial cells. *Int J Mol Med*. 2013 Feb;31(2):493-7.
33. Li D, Liu Y, Yang Y, Chen JH, Yang J, Zou LY, Tian ZQ, Lv J, Xia PY. Looped Host Defense Peptide CLP-19 Binds to Microtubules and Inhibits Surface Expression of TLR4 on Mouse Macrophages. *J Immunol*. 2013 Jun 15;190(12):6083-92.
34. Li L, Ye XP, Lu AZ, Zhou SQ, Liu H, Liu ZJ, Jiang S, Xu SY. Hyperglycemia magnifies bupivacaine-induced cell apoptosis triggered by mitochondria dysfunction and endoplasmic reticulum stress. *J Neurosci Res*. 2013 Jun;91(6):786-98.
35. Zhao L, Chen J, Cheng L, Wang X, Du J, Wang F, Peng Z. Effects of *Enterococcus faecalis* lipoteichoic acid on receptor activator of nuclear factor- κ B ligand and osteoprotegerin expression in periodontal ligament fibroblasts. *Int Endod J*. 2014 Feb;47(2):163-72.

Version 2016.12.08